



## BRAND HandyStep® S Dispenser



Ideal for routine and research applications in diagnostics, molecular biology, environmental analysis, and many more. HandyStep® S and PD-Tips work on the direct displacement principle. This offers the highest-precision dispensing of liquid media with high viscosity, high density, or high vapor pressure. Direct displacement permits contamination-free operating since no aerosols are formed. Supplied with conformity-certification, performance certificate, operating manual, shelf/rack mount, 0.1, 1 and 10mL tips.

- Serial pipetting made easy
- Offers 59 different partial volumes with different numbers of dispensing steps, depending on tip size and stroke setting number used
- Easy tip mounting
- Optimal results are obtained with PD-Tips from BRAND, also suitable for use with Encode™ tips, Repet tips, Combitips®, Combitips® plus and other compatible tips
- Robust, stable and durable
- Dispensed volume range from 2µL to 5mL
- Increased chemical resistance through innovative plastic materials

Setting	Tip Size, mL										Steps
	0.1	0.5	1	1.25	2.5	5	10	12.5	25	50	
1	2	10	20	25	50	100	200	250	500	1000	49
1.5	3	15	30	37.5	75	150	300	375	750	1500	32
2	4	20	40	50	100	200	400	500	1000	2000	24
2.5	5	25	50	62.5	125	250	500	625	1250	2500	19
3	6	30	60	75	150	300	600	750	1500	3000	15
3.5	7	35	70	87.5	175	350	700	875	1750	3500	13
4	8	40	80	100	200	400	800	1000	2000	4000	11
4.5	9	45	90	112.5	225	450	900	1125	2250	4500	10
5	10	50	100	125	250	500	1000	1250	2500	5000	9
Volume, µL											

Code	Alt Ref	Weight, g	Price
PIP9340	705110	108	£332.00

### Accessories

Code	Alt Ref	Description	Price
PIP9342	705130	Wall support	£20.80

## BRAND HandyStep® touch



Repetitive pipettes are the first choice when it comes to efficient dispensing of long series. The HandyStep® touch repetitive pipette automatically recognises the volume of the new BRAND PD-Tips II. This saves time and reduces the risk of error. Tips from many other manufacturers can also be used, providing maximum flexibility. Charging is easy and efficient: simply place the HandyStep® touch in the inductive charging stand and the battery automatically starts charging.

- Touchscreen operation with intuitive menu navigation in English, German, Spanish, French and Chinese
- HandyStep® touch allows multi-dispensing, auto-dispensing and pipetting; HandyStep® touch S includes additional features such as sequential dispensing, multi-aspiration and titration
- Automatic size detection of BRAND PD-Tips II
- Compatible with PD-Tips prior to the PD-Tips II generation and with tips from many other manufacturers
- Automatic tip ejection
- Variable volume adjustment from 1.0µL to 50mL
- Inductively rechargeable