



**TWIRL'EM®**

## TWIRL'EM®

The TWIRL'EM® provides a safe, contaminant-free and flexible container that ensures reliable test results. Its integrated secure closure system makes it a sealed and airtight solution. Proven efficient and versatile, it is used by many professionals in various industries.

**The simple and economical tool that will support you in all your sampling activities.**

## Features and benefits

- Made of FDA compliant material
- Can accommodate solid, semi-solid and liquid samples
- Maximum bag opening to facilitate insertion of sample
- Various sizes and closure combinations available depending on the type of sample
- Suitable for use in the homogenizer
- Guaranteed sterility
- RNase, DNase and pyrogen free




**Inspired by your challenges**

**LABPLAS**


VISIT [LABPLAS.COM](http://LABPLAS.COM) FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343



# TWIRL'EM® BAG INSTRUCTION

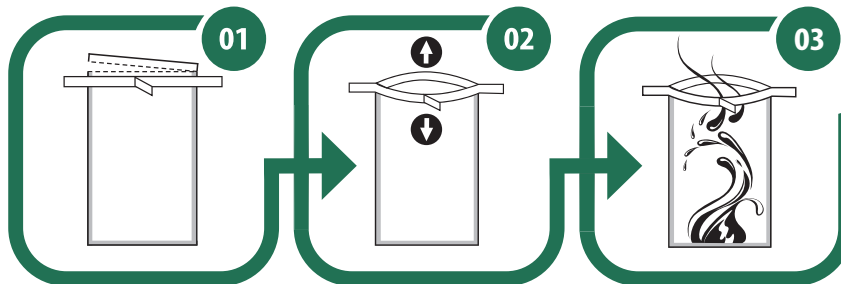

Tear off the top of the bag along the perforations

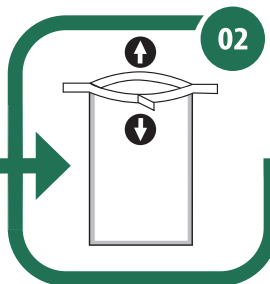


Use pull-tabs to open the bag




Aseptically insert the sample in the bag



Pull the wire-ends on each side of the bag

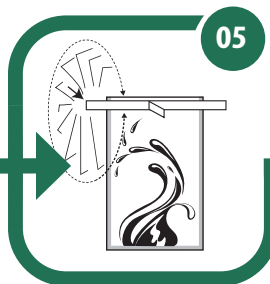

Twirl the bag around the wires, at least 3 times, to close the bag

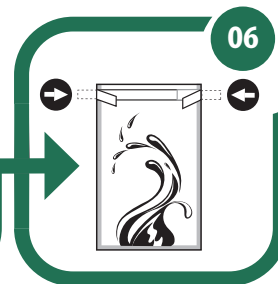

Fold both wire-ends onto the bag to provide airtight and leak-proof closure



Inspired by your challenges

LABPLAS

VISIT [LABPLAS.COM](http://LABPLAS.COM) FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343

