
Vacuum pump  
RE3022C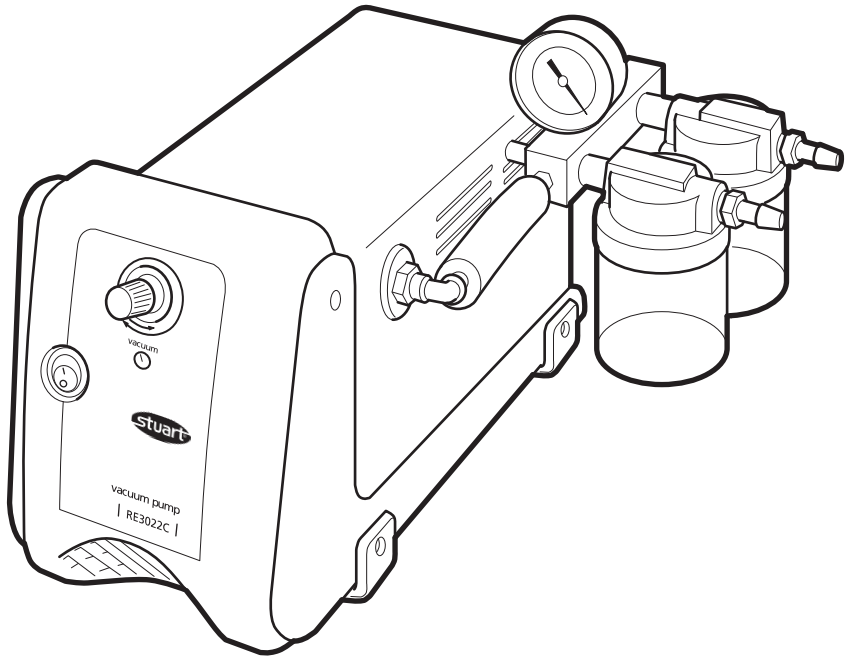

## Instruction Manual

Version 1.2

# English

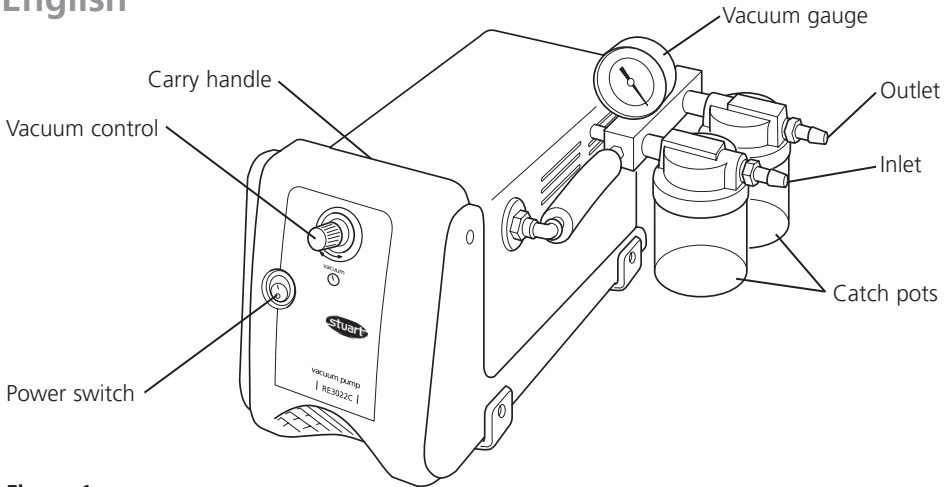
Figure 1

## VACUUM PUMP RE3022C Instructions for use

To get the best performance from the equipment and for your own safety, please read these instructions carefully before use.

If the equipment is not used in the manner described in this manual the protection provided may be impaired.

This Stuart vacuum pump is designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuations not exceeding 10% of nominal
- ❖ Overvoltage category II IEC664
- ❖ Pollution degree 2 IEC 60364-4-443
- ❖ Use with a minimum distance all round of 200mm from walls or other items

## Electrical installation

**THIS EQUIPMENT MUST BE EARTHED**

Before connection please read and understand this manual and ensure that the line supply is suitable. This pump requires a supply rated at 230V, 50Hz, single phase, ~.

Power consumption is 150W.

The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument. One has a U.K. 3 pin plug and the other has a 2 pin "Schuko" plug for connection to the mains. Choose the lead appropriate for your electrical installation and discard the other.

Should neither lead be suitable, take the lead with the UK plug and replace the plug with a suitable alternative. This involves cutting off the moulded plug, preparing the cable and connecting to the rewirable plug in accordance with its instructions.

The wires in the mains cable are coloured as follows:

- |              |           |
|--------------|-----------|
| BROWN        | - LIVE    |
| BLUE         | - NEUTRAL |
| GREEN/YELLOW | - EARTH   |

**IT IS IMPORTANT THAT THIS OPERATION IS ONLY UNDERTAKEN BY A QUALIFIED ELECTRICIAN**

The unit is fitted with an IEC socket on a flying lead at the rear of the instrument. This lead and the suitable mains lead should be connected together BEFORE connection to the mains supply.

Should the mains lead need replacement a cable of 1mm<sup>2</sup> of harmonised code H05W-F connected to an IEC 320 plug should be used.

N.B. The UK mains lead is protected by a 10A fuse mounted in the plug top.

**IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN**

## Safety advice

- ❖ If the mains supply is interrupted the units will restart with its current settings on the restoration of the electricity supply
- ❖ The pump is protected thermally. If the thermal protection has been activated the pump may restart unexpectedly when the protection resets on cooling.
- ❖ Before use ensure that the outlet port is not blocked. Before first use remove the cap fitted for protection during transit.
- ❖ When using any equipment under vacuum ensure adequate protective clothing is worn. At a minimum this should be eye protection, gloves and overalls.
- ❖ Do not operate the pump in an atmosphere containing flammable or explosive vapours.
- ❖ Do not lift or carry the pump while it is switched on or connected to the electricity supply.
- ❖ The unit should be carried using the handle located on the top (see fig 1).

## Before use

See figure 1

Before installing and using the pump for the first time please familiarise yourself with the general layout and features of the pump.

In particular ensure you have correctly identified the inlet and outlet ports. These are clearly labelled on the top of the pump heads.

The on/off switch is located at the front of the motor. Position (I) is on and position (O) is off.

There are plastic catch pots fitted to both outlet and inlet ports to collect liquid ejected from the outlet and prevent liquid ingress through the inlet. The condition of the catch pots should be checked before each use for cracks, chips or signs of corrosion. Replace pots if defects are found.

The inlet port is fitted with a vacuum gauge.

Pressure can be adjusted via the vacuum regulator located at the front of the unit.

## Installation

1. Choose a clean, dry and well ventilated area where the pump will not be exposed to moisture or flammable vapours.
2. The unit should be placed on a firm level surface.
3. The pump is fitted with 4 rubber feet which help reduce noise and creeping while the pump is in use.
4. It is recommended that flexible hose connections between the pump and the vacuum system are kept as short as possible.
5. It is recommended that a flexible hose be attached to the outlet port in order to lead any fumes which pass through the pump away from the working area, preferably into a fume hood. Make sure this hose does not become kinked and block the outlet path.
6. The tubing connectors on the inlet and outlet ports accept tubing of 9mm internal diameter.

# Operation

## Do not use this pump to pump liquids.

1. Place the pump in the required location and connect to the electricity supply.
2. Ensure that the outlet port is not blocked.

**Never block the outlet port. If the pump is run with the outlet blocked pressure will build up in the pump which can lead to the pump head bursting.**

3. Ensure that the catch pots are empty and in good condition. (see fig 1).
4. Switch the pump on and allow it to run for a few minutes in order to warm up before connection to the vacuum system.
5. Switch off and connect the pump to the vacuum system.
6. Fully open the vacuum regulator (see fig 1) by turning it anti-clockwise to the stop.

**Note:** It is necessary to completely turn the vacuum regulating valve a number of times before it is fully open. When fully open the vacuum regulator will reach a stop point and won't be able to turn any further, take care not to force the valve past this point.

7. Switch on the pump.
8. Slowly close the vacuum regulator until the required vacuum level is shown on the gauge (see fig 1).

**Do not run the pump continuously for more than 18 hours at blankoff. This will cause the pump to overheat and seize.**

9. After use release the vacuum before switching off the pump. This can be done either by venting the vacuum system or fully opening the pump's vacuum regulator.
10. Switch off the pump and disconnect from the vacuum system.

11. It is recommended that after use the pump be switched on again and run while disconnected from the vacuum system for 2 minutes. The air passing through will purge any vapours or condensate from the pump heads and extend the diaphragm life.

## Safety features

For the protection of the user and the pump there are several safety features included (See fig 1).

### 1. Inlet Catch Pot

Liquid or particulate matter entering the pump will cause damage and loss of performance. In order to prevent this a catch pot is fitted to the inlet port to trap such material before it can enter the pump head.

If the pot becomes half full it should be emptied as follows:

1. Switch off the pump and release the vacuum
2. Unscrew the catch pot
3. Empty, clean and dry and examine the catch pot and seals before screwing back into place
4. Restart the pump.

### 2. Outlet Catch Pot

If a heavy vapour load is pumped it is possible for condensation to occur inside the pump because of the compression. The pump is designed to cope with this load and will not be damaged. However the resulting liquid will be ejected from the outlet port.

The catch pot will collect this liquid preventing it being ejected into the laboratory.

If the pot becomes half full it should be emptied using the procedure outlined above for the inlet pot. The contents of the pot should be disposed of according to appropriate legislation.

## Maintenance

Periodically clean the pump with a damp cloth and a mild detergent solution

This vacuum pump is 100% oil-free. All bearings are sealed for life and permanently lubricated. Lubrication should not be attempted.

It is recommended that any servicing or repair is only undertaken by suitably qualified personnel.

Only spare parts supplied by Cole-Parmer or its agent should be used. Fitting of non-approved parts may affect the performance of the safety features of the instrument.

If in doubt, contact the Technical Department of Cole-Parmer.

Tel: +44 (0)1785 810433

The following spare parts are available from your laboratory supplier. For a comprehensive parts list please contact the Service Department quoting model and serial number.

<b>Description</b>	<b>Cat. No.</b>
Catch pot	RE3022C/1
Catch pot seal	RE3022C/2

## Warranty

Cole-Parmer warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of three (3) years. In the event of a justified claim Cole-Parmer will replace any defective component free of charge.

This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage by installation adaptation, modification, fitting of non-approved parts or repair by unauthorised personnel.

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Cole-Parmer Ltd.  
Beacon Road,  
Stone, Staffordshire,  
ST15 0SA, United Kingdom  
Tel: +44 (0)1785 810475  
Email: cpservice@coleparmer.com  
Web: www.stuart-equipment.com

# Notes

This product meets the applicable harmonized standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in its vicinity will meet these standards

and so we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, contact the manufacturer.

## EU Declaration of Conformity

<b>Product</b>	Laboratory Equipment	<b>File Number</b>	P225
<b>Manufacturer</b>	Cole-Parmer Ltd Beacon Road Stone, Staffordshire ST15 0SA United Kingdom		

This declaration of conformity is issued under the sole responsibility of the manufacturer

**Object of Declaration** Vacuum Pumps  
(reference the attached list of catalogue numbers)

The object of the declaration described above is in conformity with the relevant Union Harmonisation Legislation:

<b>Low Voltage Directive</b>	2014/35/EU
<b>EMC Directive</b>	2014/30/EU
<b>RoHS Directive</b>	2011/65/EC
<b>Machinery Directive</b>	2006/42/EC

References to the relevant harmonised standards used or references to the other technical specifications in relation to which conformity is declared:

IEC/EN 61010-1:2010	Safety requirements for electrical equipment for measurement, control and laboratory use. Part 1: General requirements.
CISPR 11:2009	Industrial, scientific and medical equipment - Radio-frequency disturbance characteristics - Limits and methods of measurement.
IEC/EN 61326-1:2013	Electrical equipment for measurement, control and laboratory use. EMC requirements. Part 1: General requirements.

Signed for and on behalf of the above manufacturer


**Additional Information** Year of CE Marking: June 2009

**Place of Issue** Stone, Staffordshire, UK

**Date of Issue** 23 January 2017

**Authorised Representative** Steve Marriott

**Title** Technical Director

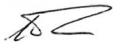**Signature** 

**Cole-Parmer Ltd.**

Beacon Road,  
Stone,

Staffordshire,  
ST15 0SA,

United Kingdom

Tel: +44 (0)1785 812121

Email: [cpinfo@coleparmer.com](mailto:cpinfo@coleparmer.com)

Web: [www.stuart-equipment.com](http://www.stuart-equipment.com)