
Cole-Parmer® HI-200 Series Hybridisation Incubator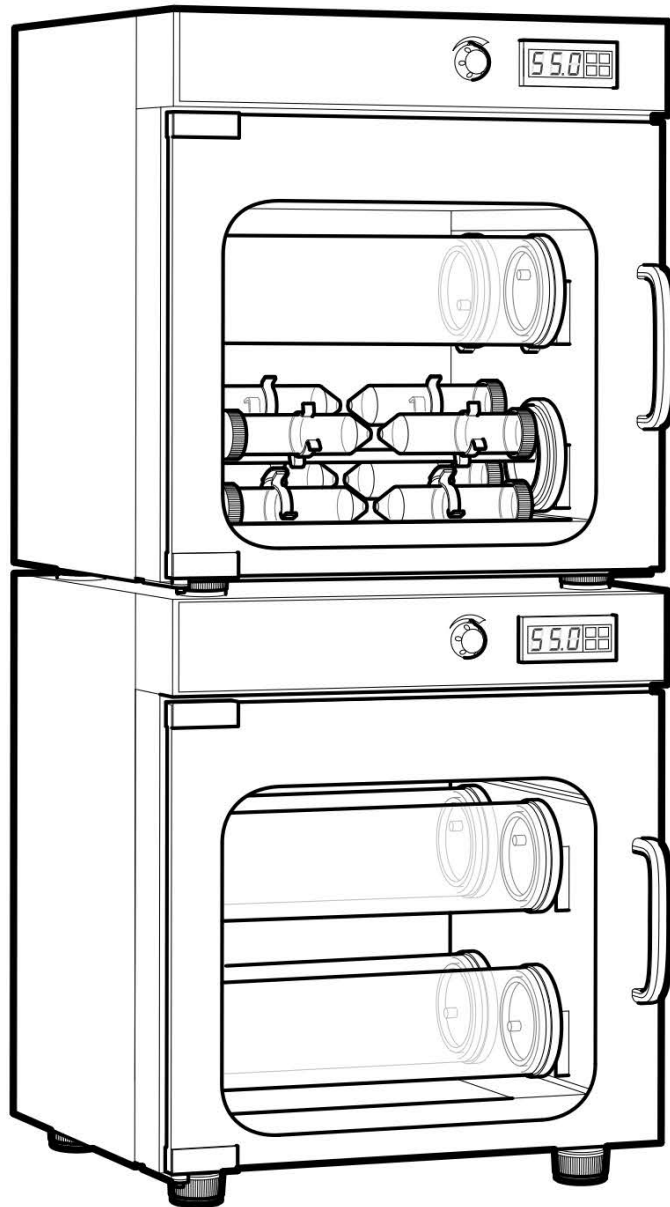

Instruction Manual
Version 8.2

Cole-Parmer®
essentials

CONTENTS

Contents

INTRODUCTION	4
BEFORE USE	5
UNPACKING	5
SAFETY AND INSTALLATION	6
WARNING	6
OPERATOR SAFETY	6
INSTALLATION	6
Replacement cable.....	7
ENVIRONMENTAL CONDITIONS	7
SÉCURITÉ ET CONSIGNES D'INSTALLATION	8
AVERTISSEMENT	8
INSTALLATION	8
Câble de rechange.....	9
CONDITIONS ENVIRONNEMENTALES	9
SICHERHEITS- UND INSTALLATIONSINFORMATIONEN	10
ACHTUNG	10
SICHERHEIT DES BEDIENPERSONALS	10
INBETRIEBNAHME	10
Ersatzkabel.....	11
UMWELTBEDINGUNGEN	11
INFORMAZIONI SULLA SICUREZZA E L'INSTALLAZIONE	12
AVVERTENZA	12
SICUREZZA DELL'OPERATORE	12
INSTALLAZIONE	12
Cavo di ricambio.....	13
CONDIZIONI AMBIENTALI	13
INFORMACIÓN DE SEGURIDAD E INSTALACIÓN	14
ADVERTENCIA	14
SEGURIDAD DEL OPERARIO	14
INSTALACIÓN	14
Cable de repuesto.....	15
CONDICIONES AMBIENTALES	15
SPECIFICATIONS	16
INSTALLATION	17
See also the sections on Safety and Installation.....	17
FITTING STACKING PLATES AND WALL MOUNTING PLATES	17
Fixing Procedure for two Hybrigenes.....	18
Fixing Procedure for three Hybrigenes.....	19
OPERATION	19
THE FRONT PANEL CONTROLS	19
SETTING THE TEMPERATURE	20
SETTING THE SPEED	20
HYBRIDISATION TUBES	21
Large (80mm diameter) tube assembly (FHB16).....	21
Small (44mm diameter) tube assembly (FHB36).....	21
Mini (4 x 32mm) tube assembly (FHB41 and FHB61).....	22
16 x 15ml tube assembly (F15ML4TH) and 8 x 50ml tube assembly (F50ML4TH).....	22
INSTALLING THE HYBRIDISATION TUBES	23

BASIC PROTOCOL FOR SOUTHERN HYBRIDISATIONS	23
Insertion of filters	23
Pre-hybridisation.....	23
Hybridisation.....	23
Washes.....	23
Notes:	23
WARNING	23
USER MAINTENANCE	24
O ring seals on the FHB16 hybridisation tubes	24
Hybridisation tubes	24
Cleaning your Hybridisation Oven.....	24
If you spill the fluid	24
ADDITIONAL INFORMATION	25
FAULT FINDING	25
FUSES	25
SAFETY OVER-TEMPERATURE CUT-OUT	25
SUPPORT CONTACTS.....	25
SPARES AND ACCESSORIES.....	26

INTRODUCTION

Please read all the information in this manual before using the unit.

The Cole-Parmer HI-200 Series Hybridisation Oven is ideal for blotting techniques in which RNA, DNA or protein are immobilised onto nylon or nitrocellulose filters. Several filters can be hybridised simultaneously in a single tube while the rotating design ensures volumes as low as 5ml can be used and recovered.

The HI-200 Series Hybridisation Oven has unique “slot-in” tubes with adjustable tube rotation speeds of 0 to 20 rpm; controllable to suit your application. The tubes and other accessories can be accommodated together for multiple applications and the unit can hold up to 16 mini or 4 large glass tubes. The temperature can be set from 10°C above ambient to 80°C and offers excellent temperature accuracy and uniformity.

For safety, the unit has a 10mm thick Perspex door to shield the immediate environment from radiation. In addition, protective casing and non-drip tube design minimises risk to the user. Included is a removable drip tray for easy cleaning should spillages occur.

The HI-200 Series Hybridisation Oven has adjustable feet for levelling on uneven surfaces. It is also possible to stack up to three ovens, saving on valuable laboratory space.

The operating instructions in this Operator’s Manual cover the following models of the Cole-Parmer HI-200 Series Hybridisation Oven:

Model number	Description	Voltage
HI-200	Hybridisation incubator	230V, 50/60Hz
HI-200-T	Hybridisation incubator, (including 4 large tubes FHB16)	230V, 50/60Hz
HI-200-120	Hybridisation incubator	120V, 50/60Hz
HI-200-T-120	Hybridisation incubator, (including 4 large tubes FHB16)	120V, 50/60Hz

BEFORE USE

Before using the HI-200 Series Hybridisation Oven please make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Antylia Scientific Ltd. or your supplier will be happy to assist you.

UNPACKING

When unpacking the unit please ensure that the following have been removed from the packaging:

- HI-200 Series Hybridisation Oven including drip tray
- Mains cables (UK and EU or US)
- Tube of silicon grease
- Hybridisation tubes x 4 (if included)
- Operator's manual
- Decontamination certificate

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Antylia Scientific Ltd. accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

SAFETY AND INSTALLATION

Please read all the information in this manual before using the HI-200 Series Hybridisation Oven.

WARNING

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Cole-Parmer have taken great care in the design of these units to protect operators from hazards, but operators should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS
- DO NOT use combustible substances near hot objects
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases.
- DO NOT place any liquid directly into the instrument.

OPERATOR SAFETY

All operators of Cole-Parmer equipment must have available the relevant literature needed to ensure their safety. It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Antylia Scientific Ltd. the protection provided by the equipment to the operator may be impaired.

All Cole-Parmer instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out.

If a safety problem should be encountered, switch off the unit at the mains socket and remove the plug from the electricity supply.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Cole-Parmer instruments are supplied with a power cable; this may be integral or plug-in.

2.  Before connecting the instrument to the mains electricity supply, check the voltage against the rating plate (located on the back of the unit). **Please note that the unit must be earthed to ensure proper electrical safety.** Connect the mains cable to a suitable plug according to the table below.

Connections	230V	120V
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

UK ONLY: The fused plug supplied with the mains cable is fitted with a 5 amp fuse to protect the instrument and the operator.

Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%.

3. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
4. Plug the mains cable into the socket on the back of the instrument.
5. Switch on the instrument. Symbols on or near the power switch of the unit have the following meanings:

- I Mains Switch On
- O Mains Switch Off

Replacement cable

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

ENVIRONMENTAL CONDITIONS

The HI-200 Series Hybridisation Oven is designed operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Relative humidity up to 80%
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2
- Use with a minimum distance all-around of 200mm from walls or other items

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of between 18°C and 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

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SPECIFICATIONS

Tube capacity	4 (large and small in any combination) 16 mini tubes (when using the FHB61 mini tube holder) 64 x 15ml tubes (in holder) 32 x 50ml tubes (in holder)
Maximum temperature*	80°C
Minimum temperature	10°C above ambient
Temperature selection	Digital
Adjustable rotation speed	0; 5 to 20 rpm
Adjustable rocking platform (optional)	30 opm recommended
Temperature display	LED
Temperature set point resolution	0.1°C
Temperature sensor	PRT
Temperature stability in the chamber	1.0°C
Temperature stability in the tubes	0.1°C
Temperature uniformity in the chamber#	±1.5°C
Temperature uniformity in the tubes#	±0.5°C
Temperature accuracy	±0.3°C
External dimensions (W x D x H)	355 x 383 x 432mm
Electrical supply	230V, 50/60Hz 120V, 50/60Hz
Power	750W
Net weight	15kg
Shipping weight	21kg

Note: The HI-200 Series Hybridisation Ovens include a pre-set over-temperature cut-out.

#Temperature measured in the centre of each tube.


*Large tubes are fitted with seals suitable for use up to 70°C. For higher temperatures, seals part number 6101488 are required.

For the hybridisation tubes see the list under Spares and Accessories

INSTALLATION

See also the sections on Safety and Installation.

1. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet and outlet vents are free from obstruction.
2. Place some water in a hybridisation tube and fit the tube in the HI-200 Series Hybridisation Oven. Adjust the feet on the base of the unit by unscrewing them anticlockwise slightly until the unit is level. Check the level by placing the tube in different positions in the oven.
3. Plug the mains cable into the power socket on the rear of the unit.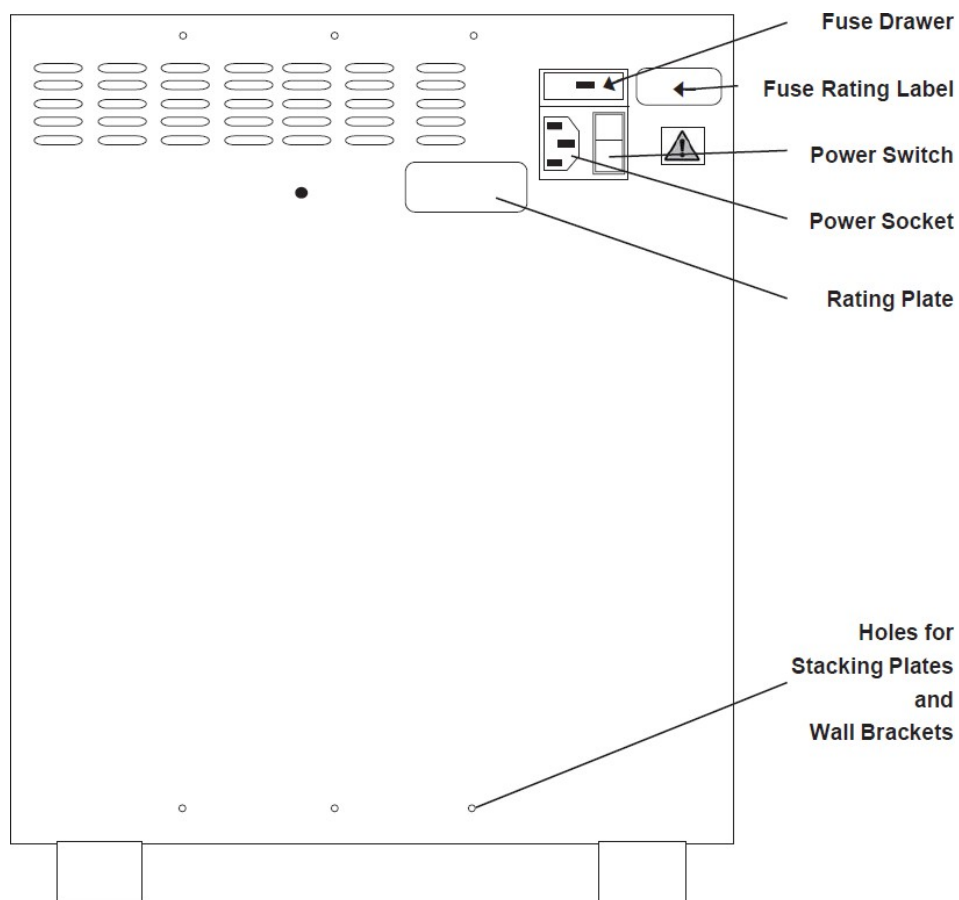

Rear view of the Hybridisation Oven

FITTING STACKING PLATES AND WALL MOUNTING PLATES

A trained technician should carry out this work only after reading the full instructions.

There are two kits available for stacking and wall fixing either two or three Hybridisation units:

FHB4DSTK: Stacking and wall mounting kit for two units


Description	Part code	Quantity
Wall Bracket	6105214	1
M4 x 8 Pozi Pan Screw	6008652	9
M4 Crinkle Washers	6008639	9
Thumbscrew	6009552	1
Wall Plate	6105213	1
Stacking Plate	6105212	1

FHB4DKIT: Stacking and wall mounting kit for three units

Description	Part code	Quantity
Wall Bracket	6105214	1
M4 x 8 Pozi Pan Screw	6008652	15
M4 Crinkle Washers	6008639	15
Thumbscrew	6009552	1
Wall Plate	6105213	1
Stacking Plate	6105212	2

Fixing Procedure for two Hybrigenes


1. Ensure the units are turned off and remove the power to them.
2. Place one unit on the top of the other.
3. The stacking plate has three holes and three slots. It has three slots to allow for the adjustment of the feet for levelling the units.
4. Use three of the screws to attach the plate to the bottom of the top unit.
5. Use another three screws to attach the plate to the bottom hybridiser.
6. Ensure that the units are level and the screws are tight.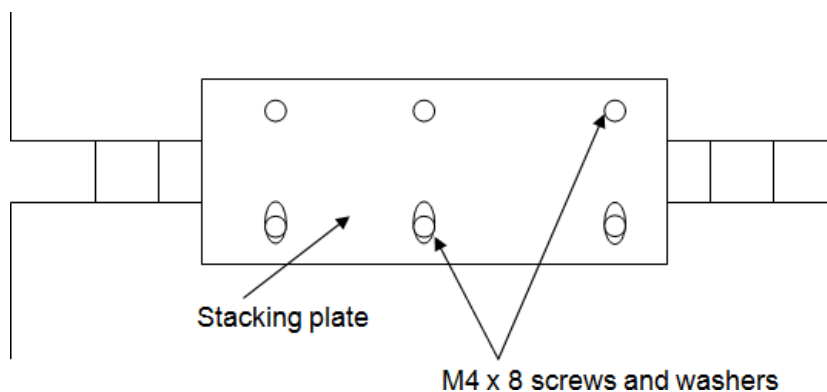

Fixing the stacking plate

7. For safety it is recommended that the top unit is fixed to a wall to prevent toppling. Use the final three screws to attach the wall bracket to the top of the top unit. Ensure that the screws are tight.

8. Put the hybridisers in the position you want to use them.
9. Position the wall plate and mark the position of the holes.
10. Use the appropriate screws and fixings to attach the wall plate to the wall; ensure that the screws are tight.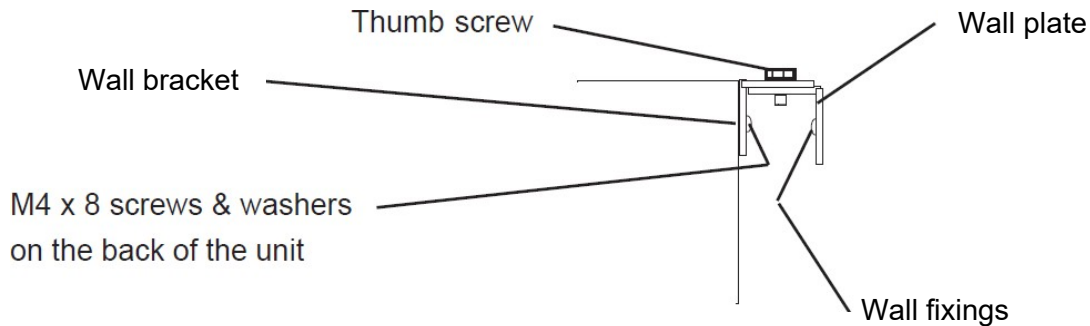

Hybridisation Oven with wall bracket attached

11. Re-position the hybridisers placing the wall bracket over the wall plate.
12. Use the thumb-screw to attach the wall bracket to the wall plate.

Fixing Procedure for three Hybrigenes

If you have three units, repeat the above process attaching a stacking plate between each pair of units.


OPERATION

SWITCHING ON

Once your Hybridisation Oven is connected to a power supply, switch it on using the power switch next to the power inlet socket on the rear panel.


THE FRONT PANEL CONTROLS

The front panel controls consist of a speed control knob for adjusting the tube rotation speed and a temperature controller.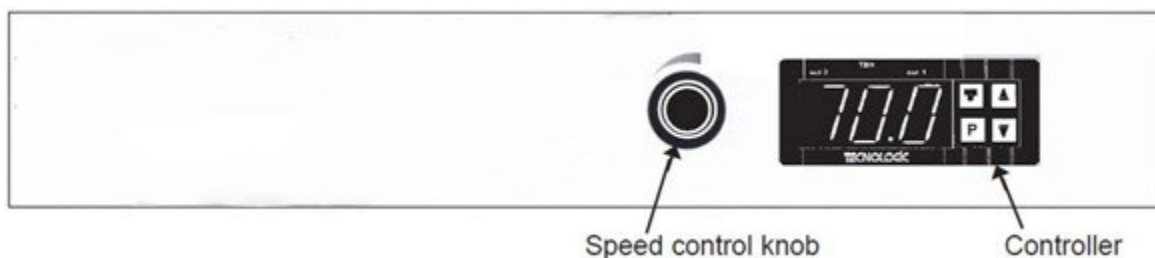

The front panel controls

SETTING THE TEMPERATURE

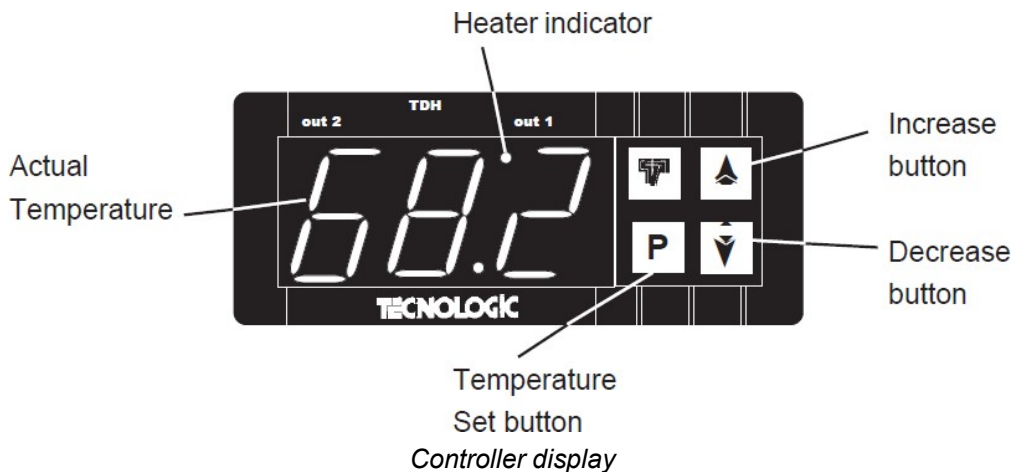1. Press the 'P' button on the controller and the set temperature will be displayed.
2. To increase the set temperature, press the Increase button.
3. To decrease the set temperature, press the Decrease button.
4. Press P and the unit will now go to the set temperature.
5. If no key is pressed for 5 seconds the display will return to show the actual temperature and the unit will go to the new set temperature.
6. The Heater indicator will be lit while the heater is on.
7. The set temperature can be checked at any time by pressing the 'P' button.
8. Press the Increase and Decrease buttons together to toggle between °F and °C.

Note: When the unit is tuned off, on subsequent power-up the units will revert to °C.

SETTING THE SPEED

Speed control knob


1. When turned fully anticlockwise, the speed control knob switches the rotation of the tubes off.
2. To increase the rotation speed, turn the knob clockwise until you reach the required speed.
3. Operation of the speed control knob increases the speed clockwise and decreases the speed anticlockwise in steps of approximately 4 rpm.

HYBRIDISATION TUBES

The hybridisation tubes are manufactured from durable thick borosilicate glass with a non-drip design. The tubes must be checked regularly and replaced if any damage occurs. Any damage to the ends of the tubes can cause them to break or leak when the end caps are inserted.

Large (80mm diameter) tube assembly (FHB16)


Remove the tube end cap which has the two holes in it. Place the filters in the tube and replace the end cap. A small quantity of Cole-Parmer O-ring grease (part code 6101351, supplied with the Hybridisation Oven) **must** be smeared on the seal to facilitate the insertion of the end caps. Remove the spindle cap and add the fluids appropriate to the protocol through the hole in the spindle using a syringe. Replace the spindle cap and press it onto the spindle so that it clicks into place. Note: the tube will not rotate properly if the cap is not fitted securely.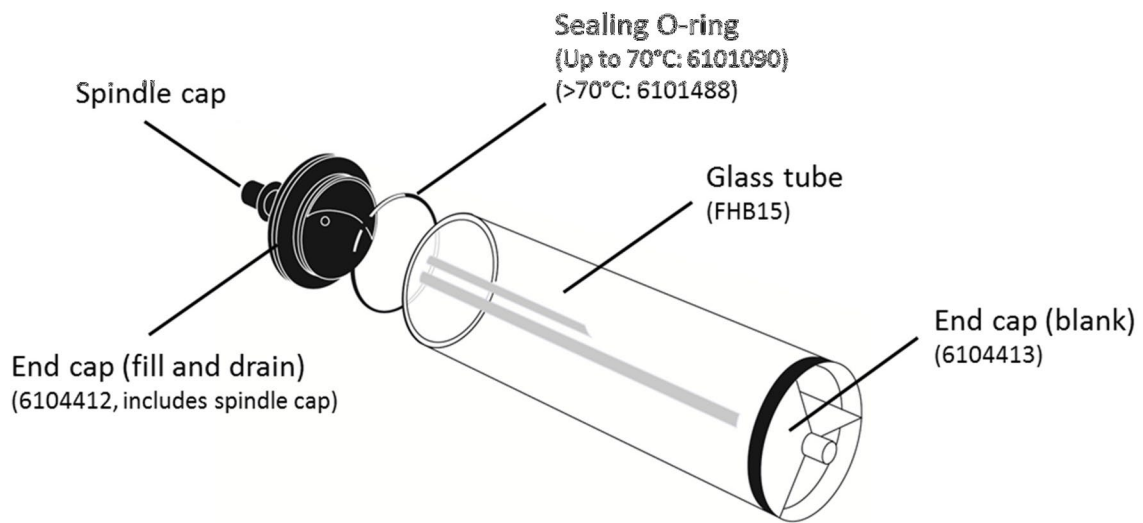

FHB16 assembly

Small (44mm diameter) tube assembly (FHB36)


Unscrew the tube end cap. Place the filters in the tube and replace the end cap. Remove the spindle cap and add the fluids appropriate to the protocol through the hole in the spindle using a syringe. Replace the spindle cap and press it onto the spindle so that it clicks into place.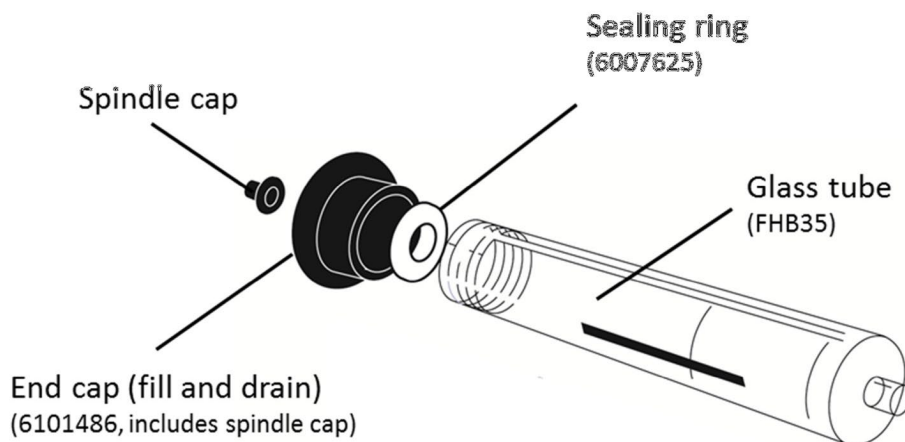

FHB36 assembly

Mini (4 x 32mm) tube assembly (FHB41 and FHB61)


The mini tubes must be used in the FHB61 multi-tube holder. Unscrew the end cap of the tube. Place the filters and the liquids appropriate to the protocol in the tube and replace the end cap. Fix the tube into a multi-tube holder, using the springs, and ensure that there are four tubes in the holder. The end caps must be as shown, on alternate ends. The multi-tube holder will not rotate properly if there are not four mini tubes in it.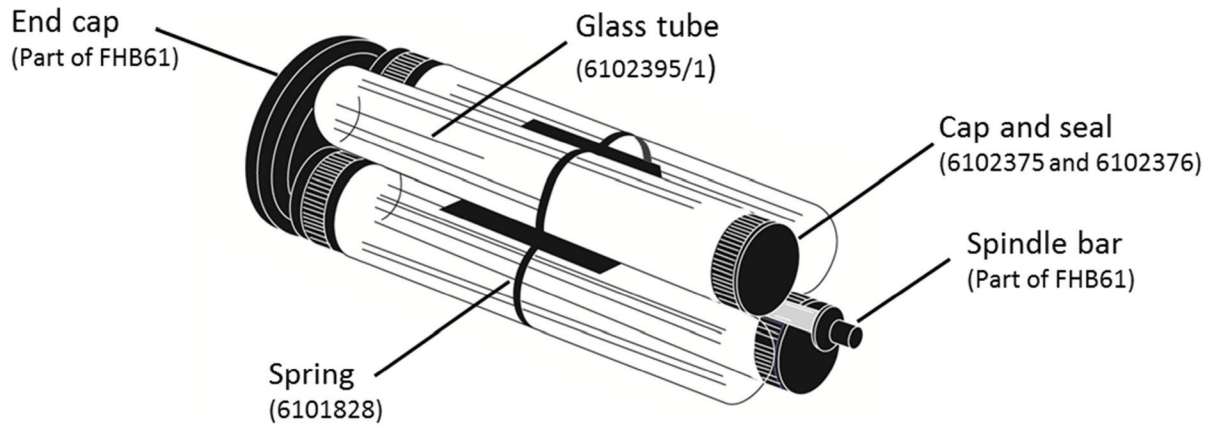

4 x FHB41 and FHB61 assembly

16 x 15ml tube assembly (F15ML4TH) and 8 x 50ml tube assembly (F50ML4TH)

These tube holders are designed to hold 15ml and 50ml Falcon-type tubes. Unscrew the end cap from the tube. Place the filters and the liquids appropriate to the protocol in the tube and replace the end cap. Assemble the tubes into the multi-tube holder using the spring clips to hold them in place. The multi-tube holder will not rotate properly if it is not balanced properly. The multi-tube holder should be filled completely or at least balanced even if all the tubes are not being used.

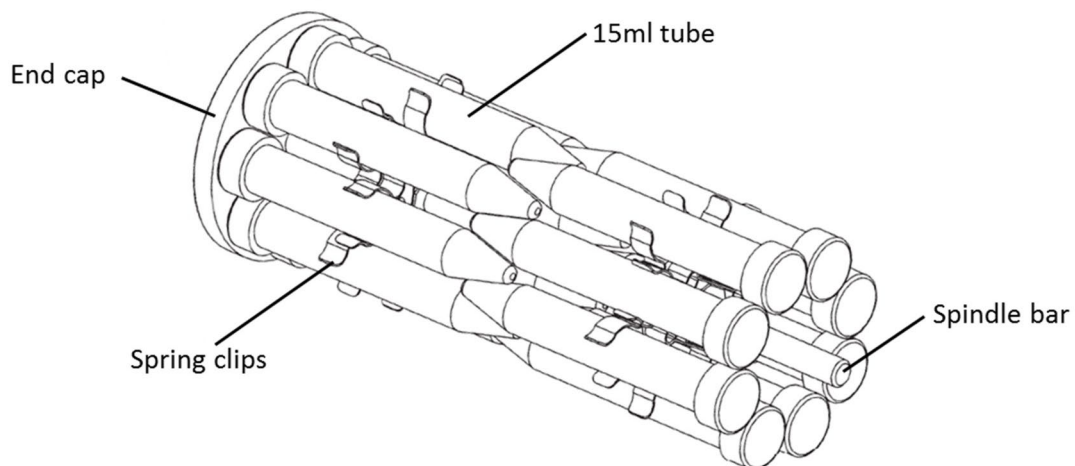F15ML4TH assembly

INSTALLING THE HYBRIDISATION TUBES

1. Pre-heat the oven to the required temperature.
2. Insert the tubes or tube holder into the Hybridisation Oven ensuring that they rotate freely and the end caps engage on the drive rollers on the right hand side of the chamber.
3. Close the door.
4. Turn the speed control knob until you reach the required speed.

BASIC PROTOCOL FOR SOUTHERN HYBRIDISATIONS

Insertion of filters

Place a wet or moist filter, DNA side facing inwards, into a hybridisation tube and roll the tube, allowing the filter to adhere to the glass. Large trapped air bubbles should be smoothed out with a glass pipette; smaller air bubbles are acceptable and do not interfere with hybridisation so long as the surface of the filter is covered with hybridisation fluid.

Several filters may be placed in one tube; there is a slight loss of signal, but no increase in background in the lower filters. To improve the signal we recommend the use of separation membranes between filters, part number FMEM2020.

Pre-hybridisation

Add 5 or 10ml of hybridisation fluid, according to the number of filters and the size of tube, so that the filters are covered. Allow to pre-hybridise for the required period.

Hybridisation

Drain off the pre-hybridisation fluid by removing the spindle cap from the fill and drain hole in the blue end cap and inverting the tube. Add the hybridisation liquid together with the radioactive probe using a syringe and needle through the fill and drain hole and replace the spindle cap. The minimum volume for a single filter is 5ml (when using large tubes). Allow to hybridise for the required period.

Washes

Drain off the hybridisation fluid through the fill and drain hole, collecting the probe in a suitable container for re-use or disposal. Washing depends on the stringency required and usually involves a series of washes with SSC buffer and SDS of decreasing concentration until background radiation has been removed.

Notes:

- All radioactive waste can be collected for disposal; none need be disposed of down the sink.
- Radioactive probes can be collected and reused several times.
- As the protocols for this unit usually involve the use of radioactive materials the appropriate manufacturer's safety precautions must be taken into consideration.

WARNING

THE UNIT WORKS UP TO 80°C. This temperature can cause burns. Antylia Scientific Ltd. takes great efforts to protect operators from hazards, but the following points are brought to your attention:

1. Use care and wear protective gloves to remove hybridisation tubes from the cabinet if the unit is at high temperature. If the protocol allows, it is better to switch off the unit and open the door to allow the unit to cool to below 40°C before attempting to remove the tubes.
2. The coloured end cap of the 80mm large tube and the 44mm small tube has two holes. Take care not to spill radioactive probe through these holes. One hole allows the air to expand when the tube is heated. The other is the fill and drain hole which is revealed when the spindle cap is removed.

DO NOT TIP THE TUBE SO THAT THE COLOURED END IS DOWN AS THIS WILL CAUSE A SPILLAGE.

3. The unit has been designed to conform to international safety specifications and is fitted with an over-temperature cut-out. If, for any reason, the temperature rises above a safe level for the instrument, the over temperature protection cuts out the heater, but the tubes continue to rotate. Check the reason for the cut out and, if there is no obvious cause, turn the unit off, allow it to cool then turn it on again. The cut-out should reset itself. If it does not re-set then the unit needs repair.
4. If the Controller displays 'Err', the thermistor has broken or become disconnected and the unit will cool down automatically.

USER MAINTENANCE

O ring seals on the FHB16 hybridisation tubes

The O ring seals on the large 80mm diameter hybridisation tubes must be checked regularly and replaced before a leak occurs. Each tube is provided with 6 spare O rings and further O rings can be obtained from Antylia Scientific Ltd. or your Cole-Parmer dealer.

Hybridisation tubes

The hybridisation tubes must be checked regularly and replaced if any damage occurs. Any damage to the ends of the tubes can cause them to leak or break when the end caps are inserted. Further tubes can be obtained from Antylia Scientific Ltd. or your Cole-Parmer dealer.

Cleaning your Hybridisation Oven

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool.

You can clean the case of the Hybridisation Oven with a cloth dampened in soapy water. No part of the unit should be immersed in solvents. Do not use acetone or abrasive cleaners.

DO NOT USE ETHANOL OR METHYLATED SPIRITS ON THE ACRYLIC DOOR; these may make the acrylic go cloudy.

Before using any cleaning or decontamination method except those recommended in this manual, the responsible body should check with Antylia Scientific Ltd. that the proposed method will not damage the equipment.

If you spill the fluid

Because of the nature of operation of the Hybridisation Oven, the complete inside must be cleaned if a spillage does occur. This is to remove any airborne particles which may have been blown around the system and to guarantee full operator safety.

Switch the Hybridisation Oven off and disconnect it from the power supply. Follow this sequence of operations when the Hybridisation Oven has been allowed to cool below at least 50°C:

- Remove all glassware from the Hybridisation Oven.
- Remove the plastic drip tray.
- Clean all surfaces of the inner chamber using a suitable cleaning and decontaminating solution. Be careful of the drive shafts.

Cleaning and decontamination materials recommended by Antylia Scientific include:

- Lipsol®
- Decon® 90
- Neutracon®
- Other mild detergent

ADDITIONAL INFORMATION

Note that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal mains voltages. There are no user serviceable parts within this equipment.

FAULT FINDING

Should you have any problems with your HI-200 Series Hybridisation Oven which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Contact the Antylia Scientific Ltd. Service Department (details given below) to obtain information on how to return the unit. Fill in the Decontamination Certificate provided with the unit and include the serial number and details of the fault observed. Please remember to return the unit in its original packing. Antylia Scientific Ltd. accepts no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit.

FUSES

If the controller fails to light when power is supplied to the unit the fuses may have blown. Check that there is no external cause such as a faulty plug; check both fuses and replace any faulty fuse with one of the correct value. Fuse values are given on the label next to the mains input socket.

SAFETY OVER-TEMPERATURE CUT-OUT

In the event of no heater power, check the mains plug and lead, then reset the cut-out control by switching the unit off, allow to cool then turn it on again. The cut-out should reset itself. If it does not re-set then the unit needs repair.

Repeated operation of the cut-out indicates a serious fault: you may need to return your unit to your supplier for repair.

SUPPORT CONTACTS

If you require further technical or application assistance please contact Antylia Scientific Ltd at:

E-mail: cptechsupport@antylia.com.

Phone: +44 (0)1785 810433

For servicing information please contact:

E-mail: cpservice@antylia.com. Phone: +44 (0)1785 810475

Antylia Scientific Ltd.
Service Department
Beacon Road
Stone
Staffordshire
ST15 0SA

Please note that neither Antylia Scientific Ltd. nor any of their agents can accept the return of ANY GLASSWARE from these units.

If the unit is returned it must be thoroughly decontaminated and a Decontamination Certificate returned with the unit. A copy of the certificate is available from our website, www.coleparmer.com, or can be requested from the Service Department if the unit is being returned.

SPARES AND ACCESSORIES

The following spares and accessories may be purchased from Antylia Scientific Ltd. or their dealers. All Cole-Parmer tubes are made from durable Borosilicate glass. Contact Antylia Scientific Ltd. or your Cole-Parmer dealer for further details.

Part code	Description
FHB16	Large (80mm diameter) hybridisation tube assembly complete with sealing rings and end caps.
FHB15	Large (80mm diameter) glass tube only
6104412	End cap (fill and drain) for large tubes, includes spindle cap (no sealing ring)
6104413	End cap (blank) for large tubes (no sealing ring)
6101387	Spindle cap (for use with 6104412 and 6101486)
6101090	Sealing ring for large tubes (for 70°C operation)
6101488	Sealing ring for large tubes (for >70°C operation)
6101351	O-ring grease (100g tube)
FHB36	Small (44mm diameter) hybridization tube assembly complete with sealing ring and end cap.
FHB35	Small (44mm diameter) glass tube only
6101486	End cap (fill and drain) for small tubes, includes spindle cap (no sealing ring)
6007625	Sealing ring for small tubes
FHB41	4 x 32mm diameter mini tubes
6102395/1	Mini tube, glass tube only
6102375	Cap for use with the mini tube
6102376	Seal for use with the mini tube
FHB61	Multi-tube holder for use with FHB41 (holds 4 mini tubes)
F15ML4TH	16 x 15ml tube assembly (holds 15ml Falcon-type disposable tubes)
F50ML4TH	8 x 50ml tube assembly (holds 50ml Falcon-type disposable tubes)
FHB4/PLAT	Rocking platform
FHB4SH1	Stainless steel mesh shelf
FHTRACK	Tube rack/holder (holds 3 large and 3 small tubes)
FMEM2020	Membrane separators 20 x 20cm (pack of 5)
FHB4DSTK	Stacking and wall mounting kit for two Hybrigenes
FHB4DKIT	Stacking and wall mounting kit for three Hybrigenes
6105112	Drip tray
HH179(S)	UK mains lead with plug, 230V
HH180(S)	European mains lead with plug, 230V
FCABLEUS	US mains lead with plug, 120V

This product meets the applicable CE Directives and UKCA Legislation for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in its vicinity will meet these standards and so we cannot guarantee

that interference will not occur in practise. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advise before use, contact the manufacturer.

Declaration of Conformity is available to view online at www.coleparmer.com

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Warranty Registration

Cole-Parmer[®]
essentials

Antylia Scientific Ltd
Beacon Road,
Stone,
Staffordshire,
ST15 0SA,
United Kingdom

UK

T: +44 (0) 1480 272279
E: uk.sales@antylia.com
W: coleparmer.co.uk

Germany

T: +49 (0) 9377 92030
E: de.sales@antylia.com
W: coleparmer.de

France

T: +33 (0) 1486 37800
E: fr.sales@antylia.com
W: coleparmer.fr

Italy

T: +39 (0) 284349215
E: it.sales@antylia.com
W: coleparmer.it

India

T: +9122 61394444
E: info@coleparmer.in
W: coleparmer.in

China

T: +1 847 549 7600
E: sales@antylia.com
W: coleparmer.com

USA

T: +1 847 549 7600
E: sales@antylia.com
W: coleparmer.com

Canada

T: +514 355 6100
E: info@antylia.ca
W: coleparmer.ca

Other

T: +1 847 549 7600