Lovibond® LC 100

Quick Start Guide

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Guia rápido

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CE Declaration

Manufacturer's Name/Authorized Representative:

The Tintometer Ltd • Lovibond House • Solar Way • Solstice Park • Amesbury • UK

Phone: +44 (0)1980 664800 • Fax +44 (0)1980 625412

X-Rite, Incorporated • Siemensstraße 12b • 63263 Neu-Isenburg • Germany

Model Name: LC 100

Model No.: LC 100

Directive(s) Conformance: EMC 2004/108/EC LVD 2006/95/EC

Federal Communications Commission Notice

NOTE: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Industry Canada Compliance Statement

This Class A digital apparatus complies with Canadian ICES-003.

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Equipment Information

Use of this equipment in a manner other than that specified by The Tintometer Ltd may compromise design integrity and become unsafe.

WARNING: This instrument is not for use in explosive environments.

This document is intended to familiarize you with the Lovibond® LC 100 instrument.


Table of Contents

1. Packaging Contents	2
2. Instrument Description	3
3. Getting Started	4
Charging the Battery	4
Powering On and Off (battery operation only)	4
Language Selection	5
Instrument Calibration	5
Sample Measurement and Preview	5
4. Function Screens	6
5. Creating Standards	7
6. Colour QC	8
Measuring Samples	8
Saving a Tag with a Measurement	8
7. Outputting Reports	10
USB Report Mode	10
8. Quick Check Menu	11
9. Sample Records Menu	11
10. Appendices	12
Troubleshooting	12
Screen Messages	12

1. Packaging Contents

- LC 100 instrument
- Lanyard
- USB interface cabling
- Protective zipper carry case
- Protective pouch with clip

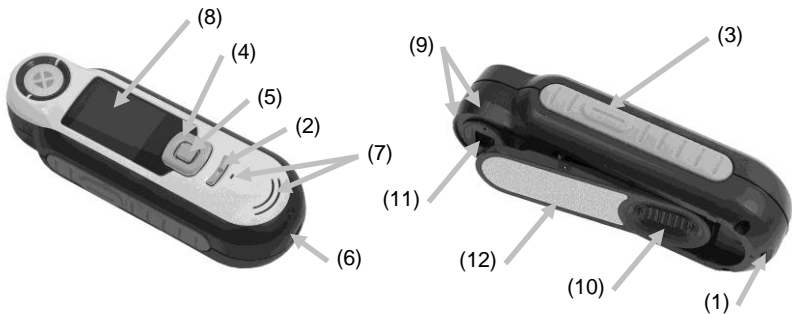2. Instrument Description

1. **Lanyard loop:** Provides a secure way to carry your device.
2. **Power on/off:** Used to turn the instrument on and off.
3. **Measure button:** Two stage button; press lightly to start preview and firmly to measure. Lightly pressing this button always returns you to the Colour QC screen.
4. **Navigation control:** Navigates the on screen menus and function bar in four directions. Pressing on the left side moves the focus to the next available item to the left. Pressing on the right side moves the focus to the next available item to the right. The top and bottom sides perform the same function, only in an up and down direction.
5. **Enter button:** Selects menu items and opens tag menu.
6. **USB connection:** For charging the battery and interface to the software.
7. **Speaker and Microphone:** For recording and playback of voice tags.
8. **Display:** Communicates color information, instrument status, and option.
9. **Targeting marks:** Allows for easy positioning on samples.
10. **White Reference slider:** Positions the calibration tile for measurement or calibration/storage.
11. **Measurement optics:** Area that illuminates and measures the color.
12. **Information label:** Provides compliance, serial number and other information.

3. Getting Started

Charging the Battery

Connect the USB cable to the LC 100 and your computer's USB port. A battery charging screen momentarily appears. For the first time connection, a language selection screen also appears. Refer to the Initial Startup section on the following page for first time operation.

Note: Some computers have low power USB ports which are typically located at the front. It is recommended that the LC 100 is **ONLY** connected to high power ports.

During charging, the instrument switches on and indicates charging with a "lightning bolt" in the battery icon. The LC 100 will not power off while plugged into the computer USB port. A full charge from empty will take 6 hours connected to a computer USB port.

Powering On and Off (battery operation only)

Turn the instrument on by pressing and releasing the **Power** button.

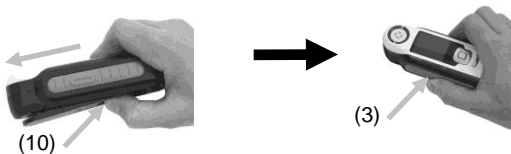
Language Selection


Select the language from the list by pressing the up and down sides of the **Navigation** control. Refer to the Introduction section for button locations.

Instrument Calibration

1. To calibrate, position reference slider (10) over optics and press **Measure** (3).

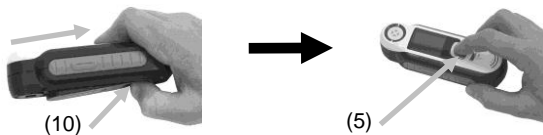
2. When finished, position reference slider (10) back to storage location and press Enter (5).


Sample Measurement and Preview

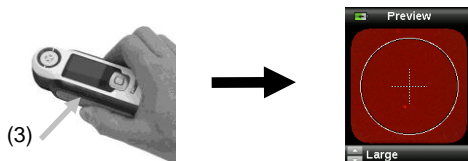1. To measure, lightly press and hold **MEASURE** button (3) and position instrument on sample.

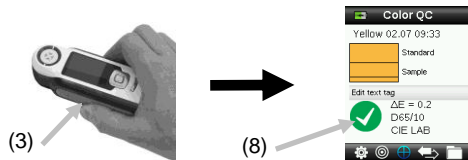
2. Press **MEASURE** firmly (3) to take measurement. View results (8).

4. Function Screens

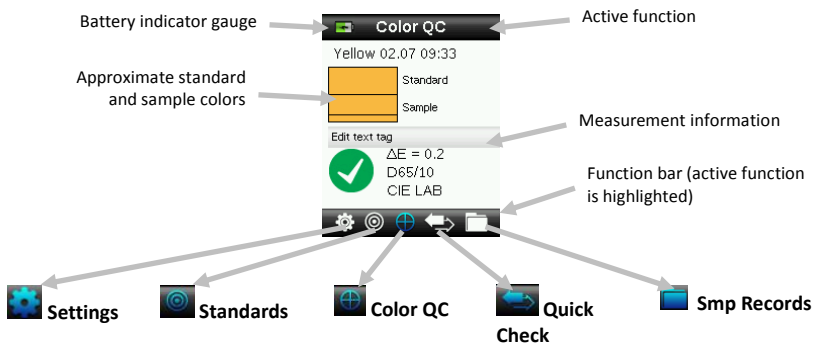Functions are selected by pressing the left or right sides of the **Navigation** control. The active function is indicated by the highlighted icon on the function bar and the title of the window.




Configure the instrument.

Select from a list or measure a new standard.


Perform Colour QC sample measurements.

Perform Colour Compare and Opacity measurements.

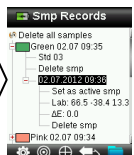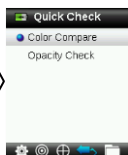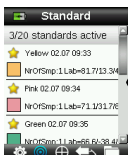
View and delete saved sample records.

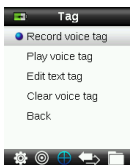
Press **Enter** (tag highlighted) Press **Navigation** control (bottom)

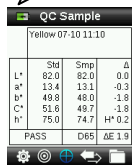Tag menu is for adding notes to a sample measurement.

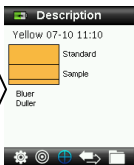Sample image, Colour plot, and Indices views are also available.


5. Creating Standards

1. Select the Standard icon in the function bar by pressing on the right or left side of the **Navigation** control.

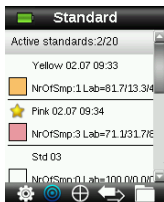2. Press the up and down sides of the **Navigation** control to select the next available standard location.

Standards active for auto selection

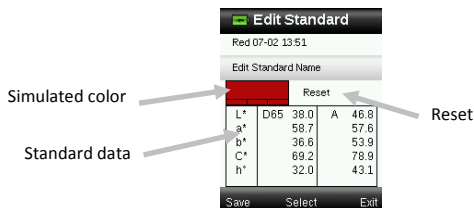
Star indicates selected standard for QC measurements

Available standard location

1. Press the left side of the **Navigation** control (Edit) to enter the Edit Standard screen. If your device has security enabled you will be prompted to enter the password “999” to unlock.
2. Preview your sample by lightly pressing then firmly to **Measure**.

Note: If you do not want the current measurement to be used, use the **Navigation** control to select Reset and press the **Enter** (Select) button.

3. Continue with additional measurements to obtain an average value if standard averaging is activated in the Settings (default setting is 3).

4. When finished, press the left side of the Navigation control (Save) to save the new standard.

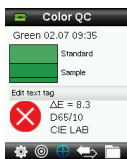
6. Colour QC

Measuring Samples

1. Select the Standard icon in the function bar by pressing on the right or left side of the **Navigation** control.

2. If the selected standards (indicated by the star) are not the standards required for sample comparisons, press the down side of the Navigation control and select the desired standard. Press the Enter button to toggle between selected (star icon) and not selected (no star icon). Continue for all standards.
3. When finished, press the right side of the Navigation control (Exit) twice to advance to the Colour QC screen.
4. Lightly press and hold the Measure button to activate the preview. The area to be measured is indicated in the preview by the circle.
5. When you have located the area you wish to measure, press the Measure button firmly.
6. Continue with additional measurements to obtain an average value if sample averaging is activated in the Settings.
7. After the measurement, a simulated colour patch of the sample appears along with the Pass/Fail indication. The sample below would indicate that the sample failed based on the tolerance set in the Setting menu.


8. Press the down side of the Navigation control to advance to additional QC Sample data screens.

Saving a Tag with a Measurement

Each measurement is automatically saved in the sample record. A voice tag or text tag can be saved with each sample measurement.


The LC 100 can store a maximum of 350 sample measurements which are stored in date & time order. When the limit of 350 stored records is reached, the oldest record is automatically deleted.

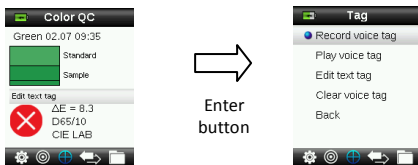1. From the Colour QC window, press the **Enter** button to display the Tag window.


2. **Voice Tag**

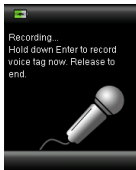To record a voice tag for the active color, select **Record voice tag** from the list and press and hold the **Enter** button down. Speak into the microphone located below the power button. When finished, release the **Enter** button.

3. **Text Tag**

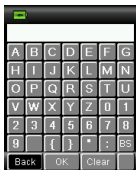To add a text tag for the current sample, select **Edit text tag** from the list and press Enter. Add your note with the virtual keyboard using the **Navigation** control and **Enter** button. When finished, highlight **OK** on the virtual keyboard and press the **Enter** Button.

4. Exit the Tag window by selecting **Back** from the list and pressing the **Enter** button.

7. Outputting Reports

USB Report Mode

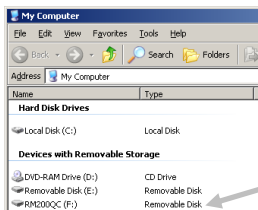
The USB report mode option is used to set the device USB mode for output of reports to computer. When the device is in this mode, it will work with your computer with similar function to a USB memory stick.

1. Plug your device into the USB port of your computer.
2. Select the USB Report Mode option from the list by pressing the up and down sides of the **Navigation** control.

3. Press the **Enter** button to switch the device in to the USB Report mode. The device will display the USB Report Mode Active window

4. The device will now be recognized on your computer as a removable disk "LC 100". If this is the first time you have plugged in your device running in the USB Report mode, the new device installation prompt will be seen and the required drivers will automatically be installed. Double click the LC 100 icon to explore the contents.

5. Two report options are provided:
 - a. QCREPORT.PDF – This provides a multipage formatted report ready for you to use directly for your quality records.
 - b. DATA.CSV – This provides data for the measured standards and samples that can be easily manipulated in Excel to a format of your preference or for more advanced data collection and analysis.

8. Quick Check Menu

The Quick Check menu contains functions which allow you to make quick comparison measurements that are not stored to the instrument. The following functions can be selected from the menu.

- Colour Compare is a quick function to measure a standard and then a sample and display results.
- Opacity Check displays the percent opacity of a drawn down sample on an opacity card.

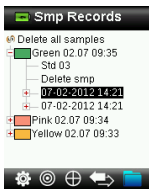
9. Sample Records Menu

The Sample Records function enables you to retrieve a stored sample from the instruments database, delete an individual sample, delete all stored samples.

1. Select the Sample Records icon in the function bar by pressing on the right side of the Navigation control.

2. Scroll through the stored records by pressing the up and down sides of the Navigation control.
3. Samples are stored under the standard that was used for measurement. Select the desired standard and press the **Enter** button to view stored sample measurements.

10. Appendices

Troubleshooting

Symptom	Action
Will not turn on	Press and hold the Power button. Charge battery / plug in to USB port.
Instrument not responding or instrument will not switch off.	1. Unplug the USB connection. 2. Hold down on/off button 10 seconds to turn off then restart.
Too much ambient light detected	Make sure you keep the instrument on the sample until the end of the measurement. Shield the sample to keep light from entering the optics during a measurement on very rough or curved samples. On transparent samples try to prevent light from entering the samples from sides or back. Place sample on a white background.
Instrument requests a "Recalibrate"	Recalibrate the instrument using the Calibrate function in the Settings menu.
"Reference slider not closed!" continually displays after calibrating more than once.	You will need to perform seven calibration procedures in a row. This will set a new reference point and correct the problem.
I cannot hear the voice recording.	Increase the volume in the Settings menu. Make sure you hold down the Enter button during recording.

Screen Messages

Screen messages can appear in the display during error conditions or for informational purposes. Some messages are cleared from the instrument screen by pressing the **Enter** button.

Messages may also be accompanied by a number to indicate a specific condition. If an error condition persists, contact technical support at +44 (0)1980 664800 or visit www.lovibondcolour.com.

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Conformité aux directives : EMC 2004/108/EC LVD 2006/95/EC

Federal Communications Commission Notice (É.-U.)

NOTE: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

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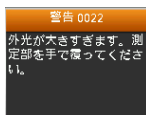
Tel: +603 3325 2286
Fax: +603 3325 2287

Tintometer China

Room 1001, China Life Tower
16 Chaoyangmenwai Avenue
Beijing, 100020
CHINA

Tel: +86 10 85251111 ext. 330
Fax: +86 10 85251001

Sales@tintometer.com

www.lovibondcolour.com

Lovibond® & Tintometer® are Registered Trade Marks of The Tintometer Limited. All translations and transliterations of LOVIBOND® & TINTOMETER® are asserted as Trade Marks of The Tintometer Limited.
Registered Office: Lovibond House, UK
Errors and Omissions Excepted. Content subject to alterations without notice.