eppendorf



Reliable Performance

Centrifuge 5804/5804 R and Centrifuge 5810/5810 R



Introducing the Eppendorf 58xx Centrifuge Family

Reliable performance

The Eppendorf approach to product development is, and always has been, about giving you more. More quality, more innovation. All of our centrifuges personalize this philosophy.

They all go beyond speed and capacity to offer you unparalleled ergonomic operation, superior temperature control and outstanding product quality. Each centrifuge is designed to operate at low noise levels for a relaxing work environment and offers soft-touch lid closure for ease of use and a low profile for easy loading and unloading of samples.

Refrigerated model centrifuges have exclusive features such as ECO shut-off and dynamic compressor control for maximum sample protection and energy savings.

New rotor generation

Eppendorf is continuously updating the centrifuge portfolio to meet changing customer needs.

Eppendorf is now introducing a new rotor generation elevating your centrifuge performance to a new level.

Enjoy unique product features that improve handling and performance. A special highlight: The new swing-bucket rotors increase the capacity of Centrifuge 5804/5804~R to $4\times250~mL$ and Centrifuge 5810/5810~R to $4\times750~mL$. Centrifuge 5810/5810~R is the most compact 3 liter centrifuge currently available.

The Benchmark in Ergonomic Operation

Low profile for easy operation

The 58xx centrifuges are designed with a particularly low height in mind. Not only does the low profile facilitate easy loading and unloading, it also allows everyone in the lab to easily reach the open centrifuge lid.



Soft-touch lid closure

There's no need to exert unnecessary high force when closing the centrifuge lid. The soft-touch lid closure allows closing the centrifuge with ease.



Eppendorf QuickLock® caps for one-hand operation

Aerosol-tight caps provide the needed safety when working with hazardous samples such as viruses, bacteria, blood or radioactive samples. The caps for the new 4×750 mL swing-bucket rotor allow for easy, one-handed operation. Its unique design allows the closed bucket to be carried securely to the biosafety cabinet.



Eppendorf QuickLock® rotor lids for ergonomic lid locking

Our aerosol-tight QuickLock rotors close with only $\frac{1}{4}$ of a turn. This not only saves you time during every day use, it also takes repetitive stress off your wrist.





Centrifuge 5804 with plate rotor A-2-DWP



Centrifuge 5804 R with rotor S-4-72 (4×250 mL)

The Centrifuge 5804/5804 R is a high speed centrifuge for medium capacity needs. It allows for molecular applications in tubes up to 250 mL and offers additional swing-bucket and fixed-angle rotors as well as deepwell plate capacity for increased versatility. Common rotors for Centrifuge 5804/5804 R include:



 4×250 mL swing-bucket rotor for blood, bacteria, cell culture centrifugation. Max. speed $3,155 \times q$ (4,200 rpm).



2-place deepwell plate rotor for plates up to 89 mm loading height. Max. speed $2,250 \times q$ (3,750 rpm).



 6×50 mL conical tube rotor with aerosol-tight QuickLock lid for tubes from 5–50 mL. Max. speed $16,639 \times g$ (11,000 rpm).



 $30 \times 1.5/2$ mL micro tube rotor with aerosol-tight lid for safe centrifugation of hazardous samples. Max. speed $20,800 \times g$ (14,000 rpm).



Eppendorf recognizes its responsibility as a global citizen. Protecting the environment is critical to the conservation of precious natural resources. The Eppendorf epGreen initiative is another example of our commitment to social responsibility. Our goal is to constantly reduce the environmental impact of our business operations and the products that we produce.

Centrifuges in general

- > Use of brushless induction motors to provide maintenance-free operation, eliminating carbon particle emissions
- > Rotors are mostly made of recyclable aluminum
- > Optimized sound design for low noise levels to benefit your work environment
- > 100 % recyclable packaging material

Centrifuge 5810/5810 R



Centrifuge 5810 with rotor S-4-104 (4 \times 750 mL)



Centrifuge 5810 R with rotor S-4-104 (4×750 mL)

The Centrifuge 5810/5810 R is a workhorse for medium to high-throughput laboratories. It combines extraordinary versatility and capacity for both tubes and plates with an extraordinary compact footprint. Common rotors for Centrifuge 5810/5810 R include:



 4×750 mL swing-bucket rotor with optional aerosol-tight Quick-Lock caps. Max. speed $3,214 \times g$ (3,900 rpm).



Aerosol-tight deepwell plate rotor for plates up to 89 mm loading height. Max. speed $3,486 \times q$ (4,500 rpm).



 6×50 mL conical tube rotor with aerosol-tight QuickLock lid for tubes from 5–50 mL. Max. speed $20,130 \times q$ (12,000 rpm).




 $48 \times 1.5/2$ mL micro tube rotor with aerosol-tight QuickLock rotor lid for safe centrifugation of hazardous samples. Max. speed $19,083 \times g$ (13,000 rpm).

Refrigerated models additionally feature

- > Dynamic compressor control for optimized cooling performance
- > CFC-free refrigerant with an ozone depletion potential (ODP) of zero
- > ECO shut-off deactivates the compressor after 8 hours of inactivity to reduce energy consumption and extend compressor life

Energy saving through new swing-bucket rotor generation*

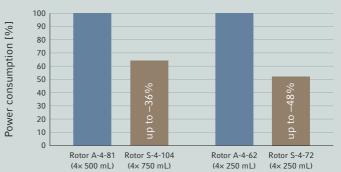

* based on internal testing with Centrifuge 5810 R, 220 V

Performance You Can Count On





> Watch video clips of Centrifuge 5810 R incl. the new rotor generation on our optimized webpage for mobile devices or www.eppendorf.com/centrifugation

Ergonomic Iid operation

> Soft-touch lid closure and low opening height for stress-free lid locking

Optimized cooling performance

> Dynamic compressor control with ECO shut-off function

Mixed loading capability

> Rotors S-4-104 and A-4-81 provide mixed loading capability with plate and tube buckets for increased versatility

Quiet operation

> All models are designed to operate at low noise levels for a relaxing work environment

Low access height (29 cm)

> Easy loading/unloading of samples

Placement flexibility

- > Compact footprint
- > Low opening height for use under shelves
- > Rolling cabinet option available for use in crowded labs

Outstanding versatility for all your applications

- > Broad rotor and adapter selection to spin all commonly used tubes, bottles, flasks, and plates
- > High centrifugation speed up to 20,913 \times g (14,100 rpm) for molecular applications in tubes from 0.2 to 750 mL
- > Aerosol-tight rotors (e.g. for potentially hazardous materials like pathogenic bacteria, viruses, blood samples or radioactive material)

The benchmark for ergonomic operation

- > Low sound level to benefit your work environment
- > Low access height (29 cm) for easy loading/unloading
- > Centrifuge lid with soft-touch lid closure

Designed to fit in every lab

- > Compact footprint saves valuable bench space
- > Low lid opening height for use under shelves
- > Automatic rotor recognition and imbalance detection for maximum operational safety

Advanced temperature management for maximum sample protection and energy efficiency

- > Temperature range from –9 °C to 40 °C
- > Maintains constant 4 °C, even at maximum speed (rotors F-34-6-38 and FA-45-6-30 maintain 4 °C at 11,000 rpm)
- > FastTemp function for quick pre-cooling
- > Standby cooling function holds temperature when centrifuge is not in use
- > ECO shut-off engages after 8 hours of non-use to reduce energy consumption and extend compressor life
- > Built-in condensation drain to eliminate water accumulation and prevent corrosion
- > Dynamic compressor control (DCC) technology for optimized cooling performance

^{*} All products which are marked with the CE-IVD symbol meet the requirements of the European Directive on IVD (98/79/EG).

Rotor Design at Its Best

Eppendorf is introducing the new generation of swing-bucket and fixed-angle rotors elevating your centrifuge performance to a new level. Enjoy unique product features that improve handling and performance.

Swing-bucket rotors

Improved energy efficiency

The new round buckets have been designed to be especially aerodynamic. The result is a more efficient design that reduces power consumption by up to 48 %.

Longer rotor lifetime

The new rotor generation of both swing-bucket and fixed-angle rotors has been designed with a particularly long lifetime in mind.

New Eppendorf QuickLock® caps for one-hand operation

- > Significantly improved handling, easy operation with one hand
- > Handle allows the closed bucket to be carried to the biosafety cabinet where it can be opened in a safe environment
- > Aerosol-tightness has been tested and verified by Public Health England, Porton Down, UK





Fixed-angle rotors

Low rotor weight

Eppendorf's fixed-angle rotors are among the lightest in the industry to improve handling and operational safety. Light-weight rotors cause less wear and tear on the motor and allow for faster acceleration and braking times.

Eppendorf QuickLock® lids for ergonomic lid locking

Our aerosol-tight QuickLock rotors close with only ¼ of a turn. This not only saves you time during every day use, it also takes repetitive stress off your wrist.

Superior temperature control

Aluminum, with its high heat conductivity, allows fast pre-cooling and accurate temperature control in refrigerated centrifuges.

Additional rotor features

- > All rotors are anodized for high chemical resistance and to prevent corrosion
- > Optimized design of rotor boreholes and adapters provide optimal support for high-speed centrifugation of tubes
- > Aluminum rotors are recyclable





Exceptional Versatility

Overview of the main rotors

Rotor image











Rotor name	S-4-104 (4 × 750 mL) aerosol-tight QuickLock caps	A-4-81 (4 × 500 mL) aerosol-tight	A-4-62 (4 × 250 mL) aerosol-tight	S-4-72 (4 × 250 mL)	A-4-44 (4 × 100 mL) aerosol-tight
5804/ 5804 R				X	X
5810/ 5810 R	Х	X	Х	X	X
Positioning and applicational range	High-capacity swing-bucket rotor for blood, bacteria, cell culture and plate centrifugation	Medium to High capacity swing-bucket rotor for blood, bacteria, cell culture and plate centrifugation	Economic swing- bucket rotor for blood, bacteria, cell culture and plate centrifugation	Economic swing- bucket rotor for blood, bacteria, cell culture centrifugation	High-speed swing- bucket rotor for platelet poor plasma, blood, bacteria and cell culture centrifugation
Aerosol-tightness	yes	yes	yes	no	yes
Max. speed (RCF)	3,214	3,220	3,220	3,234	4,500
Max. speed (rpm)	3,900	4,000	4,000	4,200	5,000
Volume range	0.2–750 mL MTP/DWP	0.2-500 mL, MTP/DWP	0.2–250 mL, MTP/DWP	0.2–250 mL	0.2-100 mL
Eppendorf Tubes® 5.0 mL and 15 mL conical	56	48	36	32	16
50 mL conical	28	20/28*1	16	16	4/8*1
13 mm blood tube	92	100	60	56	36
16 mm blood tube	80	64	48	52	28
1.5/2.0 mL micro tube	200	96	64	104	48
MTP	16	16	12		
DWP	1	1	1		

^{*1} Maximum capacity with dedicated 50 mL bucket.





Your centrifuge in safe hands with **Eppendorf Centrifuge & Rotor Service**

Quality Management Systems increasingly ask for regular inspection of centrifuges and rotors in an operational or disassembled state to ensure a safe working environment. For the instrument's peak performance and years of productive operation, an annual inspection of centrifuges and rotors by an Eppendorf certified service technician is highly recommended.



^{*2} Max. speed in Centrifuge 5810 R, lower speed applies to Centrifuge 5804/5804R/5810.

With Eppendorf service programs we make sure that temperature accuracy, rotational speed verification and operating time meet all technical specifications. Operational Qualification (OQ) is designed to safeguard your results as well as satisfy legal, governmental and internal laboratory requirements with the necessary certification documents.

Eppendorf's rotor inspection program includes cleaning, instructions for proper care and visual examination of rotors and adapters determining if the rotor is safe for further operation or needs to be replaced.

Rotor S-4-104



Tube						
Tube capacity (mL)	1.5/2.0	12 × 75 mm, 5 mL FACS	2.6-8	5.5–12	9	14
Max. tube diameter (Ø mm)	11	12	13	16	17.5	17.5
Max. vessel height (mm)	39	111	115	119	111	117
Max. vessel height for use with aerosol-tight lid (mm)*2	39	100	108	114	106	112
Adapter/bucket options for Rotor S-4-104						
Capacity per bucket/rotor	50/200	27/108	23/92	20/80	20/80	14/56
Adapter bottom shape	flat	round	round	round	round	round
Centrifugation radius (mm)	188	175	179	180	179	185
Max. RCF (g-force)	3,197	2,976	3,044	3,061	3,044	3,146
Order no. (set of 2)	5825 740.009	5825 747.003	5825 738.004	5825 736.001	5825 743.008	5825 748.000

^{*} Aerosol-tightness tested by Public Health England, Porton Down, UK.

* Exceeding the maximum tube length may reduce the loading capacity of the rotor. Please perform a manual swing-out test before starting the centrifuge run to make sure the tubes do not hit the rotor yoke.


* Please note that these conical bottles need to be used together with a conical form insert provided by the manufacturer.

* Available as a set of 2 bottles (see ordering information).

* Please note that a software upgrade will be necessary for older centrifuges when using Rotor S-4-104.

Please contact your local Eppendorf organization or dealer for further information

Order no.
5820 740.000
5820 754.001
5820 755.008
5820 743.000
5820 744.006
5820 747.005
5820 748.001
5820 708.000

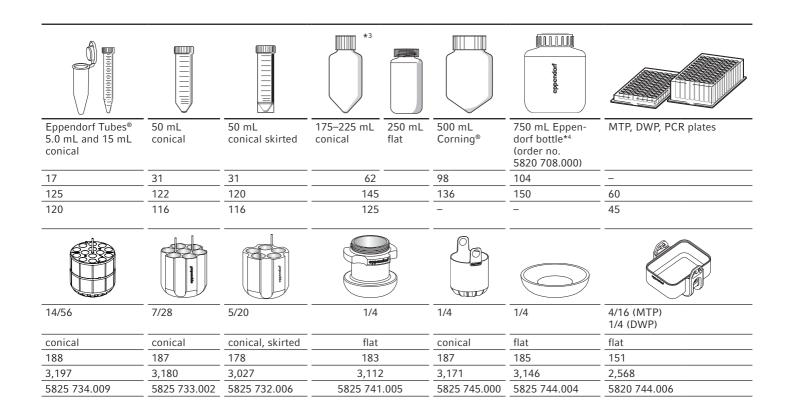


Rotor A-4-81



Tube					bre not		
Tube capacity (mL)	1.5/2	1.2-5	1.8–7	2.6–7	4–10	7–17	Eppendorf Tubes [®] 5.0 mL and 15 mL conical
Max. tube diameter (Ø mm)	11	11	13	13	16	17.5	17
Max. vessel height (mm)*2	118	118	118	118	118 (130 mm in 2 inner rows)*2	118 (130 mm in 2 inner rows)*2	122
Max. vessel height for use with aerosol-tight lid*1 (mm)	108	108	108	108	108 (130 mm in 4 center pos.)*2	108 (130 mm in 4 center pos.)*2	120
Adapter/bucket options for Rotor A-4-81							
Capacity per bucket/rotor	20/80	20/80	25/100	18/72	16/64	16/64	12/48
Adapter bottom shape	flat	flat with rubber mat	flat with rubber mat	flat with rubber mat	flat with rubber mat	flat with rubber mat	conical
Centrifugation radius (cm)	16.5	16.8	16.8	16.8	16.8	16.8	17.3
Max. RCF (g-force)	2,950	3,000	3,000	3,000	3,000	3,000	3,100
Order no. (set of 2)	5810 745.004	5810 746.000	5810 720.001	5810 747.007	5810 748.003	5810 721.008	5810 722.004
*1 Aerosol-tightness tested by Public Health Engla	nd, Porton Down, UK.						

Aerosol-tightness tested by Public Health England, Porton Down, UK.
 Exceeding the maximum tube length may reduce the loading capacity of the rotor.
 Please perform a manual swing-out test before starting the centrifuge run to make sure the tubes do not hit the rotor yoke.
 Only Centriprep® tube without locking lid.

Description	Order no.
Rotor A-4-81	
Rotor A-4-81, incl. 4 × 500 mL rectangular buckets	5810 718.007
Rotor A-4-81, incl. 4 × MTP/Flex plate buckets	5810 725.003
Rotor A-4-81 without buckets	5810 743.001
Buckets for Rotor A-4-81	
Bucket, 500 mL rectangular, set of 4	5810 730.007
Bucket, 7×50 mL conical , set of 4	5825 730.003
Bucket, MTP/Flex plates, set of 2	5810 742.005
Bucket, MTP/Flex plates, set of 4	5810 741.009
Accessories	
Aerosol-tight cap ² for 500 mL rectangular bucket, set of 2	5810 724.007
Bottle 400 mL wide-mouth, includes stabilizing sleeves, set of 2	5810 729.009
Bottle 500 mL wide-mouth, set of 2	5820 707.003

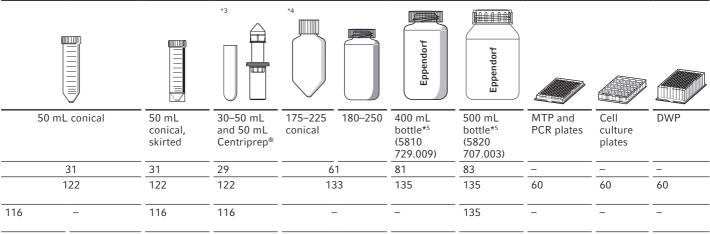













		5804 737.008			no adapter needed	See page 23 for additional ada	
5/20	7/28	5/20	1/4	1/4	1/4	4/16	2/8 1/4
conical	conical	flat with rubber mat	flat with rubber mat	flat, rubber	-	flat with rubber mat	
17.3	18.0	17.3	17.3	18.0	18.0	16.3	
3,100	3,250	3,100	3,100	3,220	3,220	2,900	
5810 723.000	5825 730.003*7	5810 739.004	5825 722.000	5810 728.002	_	5810 742.005	5810 741.009*6

^{**} Please note that these conical flasks need to be used together with a conical form insert provided by the flask manufacturer. *5 Available as a set of 2 flasks (see ordering information).

^{*6} Set of 4.
*7 50 mL conical tubes skirted need to be ordered separately. Order no. 5804 737.008 (set of 8).

Rotor A-4-62



Product features


- > Economic swing-bucket rotor with speed up to $3,220 \times g$ (4,000 rpm)
- > Rotor A-4-62 accommodates tubes and flasks from 0.2 mL to 250 mL as well as plates
- > Aerosol-tight caps*1 for safe centrifugation of hazardous samples optionally available
- > Rotor life time: max. 40,000 cycles
- > Rotor, buckets, caps and adapters are autoclavable

Tube				Na al			*3	*3
Tube capacity (mL)	1.5/2	1.2–5	1.8-7	4–10	7–17	Eppendorf Tube® 5.0 mL and 15 mL conical	7–18	18–30
Max. tube diameter (Ø mm)	11	11	13	16	17.5	17	20	26
Max. vessel height (mm)*2	122	123	121	121 (130 mm in 2 center rows)*2	118 (130 mm in 2 center rows)*2	127	126	119
Max. vessel height for use with aerosol-tight lid*1 (mm)	118	115	118	116 (130 mm in 2 inner pos.)	114 (130 mm in 2 inner pos.)	121	119	116
Adapter/bucket options for Rotor A-4-62								
Capacity per bucket/rotor	16/64	25/100	15/60	12/48	12/48	9/36	8/32	4/16
Adapter bottom shape	flat	flat with rubber mat	flat with rubber mat	flat with rubber mat	flat with rubber mat	conical	flat with rubber mat	flat with rubber mat
Centrifugation radius (cm)	17.1	17.3	17.3	17.3	17.3	17.8	17.3	17.3
Max. RCF (g-force)	3,000	3,050	3,050	3,050	3,050	3,150	3,050	3,050
Order no. (set of 2)	5810 751.004	5810 750.008	5810 752.000	5810 753.007	5810 754.003	5810 755.000	5810 756.006	5810 757.002

^{*1} Aerosol-tightness tested by the Centre for Public Health England, Porton Down, UK.
*2 Exceeding the maximum tube length may reduce the loading capacity of the rotor. Please perform a manual swing-out test before starting the centrifuge run to make sure the tubes do not hit the rotor yoke.
*3 For high g-forces we recommend using centrifuge tubes of Duran® glass.

Olu	ermu	IIIIU	rmation

Description	Order no.
Rotor A-4-62	
Rotor A-4-62, incl. 4 × 250 mL rectangular buckets	5810 709.008
Rotor A-4-62, incl. 4 × MTP plate buckets	5810 711.002
Buckets for Rotor A-4-62	
Bucket, 250 mL rectangular, set of 4	5810 716.004
Bucket, MTP plates, set of 4	5810 702.003
Accessories	
Aerosol-tight cap*1 for 250 mL rectangular bucket, set of 2	5810 710.006

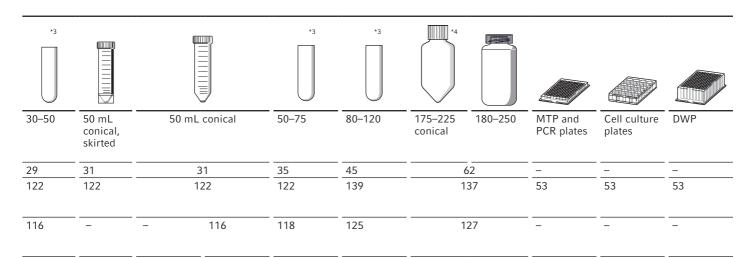















See page 23 for additional adapters

3004 /37.000								
4/16	4/16	3/12	2/8	1/4	1/4	4/16	2/8	1/4
flat with rubber mat	flat w/conical form insert	conical	flat with rubber mat	flat with rubber mat	flat with rubber mat	flat with rubber mat		r mat
17.3	17.3	17.8	17.3	17.3	18.0		15.4	
3,050	3,050	3,150	3,050	3,050	3,200	2,750		
5810 759.005	5810 763.002	5810 758.009	5810 760.003	5810 761.000	5810 770.009	5810 702.003*6		3*6

^{**} Please note that these conical flasks need to be used together with a conical form insert provided by the manufacturer.

*5 Form inserts for 50 mL conical tubes skirted need to be ordered separately. Order no. 5804 737.008 (set of 8).

*6 Set of 4.

Rotor S-4-72




Product features

- > Medium capacity swing-bucket rotor with speed up to $3,234 \times g$ (4,200 rpm)
- > Broad range of adapters available to handle all common tube formats from 0.2 mL to 250 mL
- > Aerodynamic round bucket design for up to 48 % energy savings compared to Rotor A-4-62
- > Rotor life time: max. 60,000 cycles
- > Buckets and adapters are autoclavable

Tube					
Tube capacity (mL)	1.5/2.0	2.6-8 mL	5.5–12	9	Eppendorf Tubes® 5.0 mL
Max. tube diameter (Ø mm)	11	13	16	17.5	17
Max. vessel height (mm)*1	43	115	112	113	60
Max. vessel height for use with aerosol-tight lid*1 (mm)	_	_	_	_	_
Adapter options for Rotor S-4-72		parameter		Promoted	
Capacity per bucket/rotor	26/104	14/56	13/52	12/48	8/32
Adapter bottom shape	flat	round	round	round	conical
Centrifugation radius (mm)	159	159	157	158	163
Max. RCF (g-force)	3,136	3,136	3,096	3,116	3,215
Order no. (set of 2)	5804 794.001	5804 789.008	5804 791.002	5804 792.009	5804 793.005

^{*1} Exceeding the maximum tube length may reduce the loading capacity of the rotor.
*2 Please note that these conical flasks need to be used together with a conical form insert provided by the flask manufacturer.
*3 Please note that a software upgrade will be necessary for older centrifuges when using Rotor S-4-72.
Please contact your local Eppendorf organization or dealer for further information

Description	Order no.
Rotor S-4-72*3, incl. 4 × 250 mL round buckets	5804 746.007

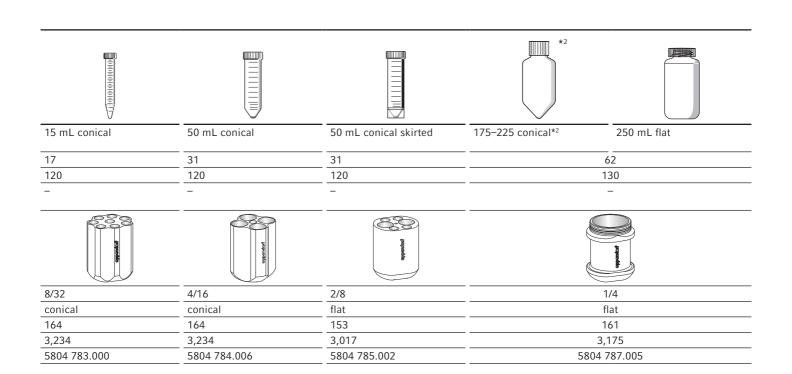

Rotor A-4-44




Product features

- > High-speed swing-bucket rotor for applications up to $4,500 \times g$ (5,000 rpm)
- > Rotor A-4-44 accommodates tubes from 0.2 mL to 100 mL
- > Aerosol-tight caps*1 for safe centrifugation of hazardous samples optionally available
- > Rotor life time: max. 34,000 cycles
- > Rotor, buckets, caps and adapters are autoclavable

Tube				lea ark			
Tube capacity (mL)	1.5/2	1.2–5	1.8–7	4–10	7–17		orf Tubes® 15 mL conical
Max. tube diameter (Ø mm)	_11	11	13	16	17.5	1	17
Max. vessel height (mm)*2	111	105	108	108	110	1	21
Max. vessel height for use with aerosol-tight lid*1 (mm)	105	102	106	106	106	_	121
Adapter/bucket options for Rotor A-4-44	0 0						
Capacity per bucket/rotor	12/48	14/56	9/36	7/28	6/24	4/16	2/8
Adapter bottom shape	flat	flat with rubber mat	flat with rubber mat	flat with rubber mat	flat with rubber mat	conical	conical
Centrifugation radius (cm)	14.8	15.0	15.0	15.0	15.0	15.5	15.7
Max. RCF (g-force)	4,100	4,200	4,200	4,200	4,200	4,300	4,400
Order no. (set of 2)	5804 751.000	5804 750.004	5804 752.007	5804 753.003	5804 754.000	5804 755.006	5804 717.007
*1Aerosol-tightness tested by Public Health England *2 Exceeding the maximum tube length may reduce Please perform a manual swing-out test before sta sure the tubes do not hit the rotor yoke.	the loading capacity of the			recommend using centrifug form inserts: 5804 728.009. 140 grams.	e tubes of Duran® glass.		

Description	Order no.
Rotor A-4-44	
Rotor A-4-44, incl. 4 × 100 mL rectangular buckets	5804 709.004
Rotor A-4-44, incl. four 2 × 50 mL buckets	5804 719.000
Buckets for Rotor A-4-44	
Bucket, 100 mL rectangular, set of 4	5804 741.005
Bucket, 2×50 mL, set of 4	5804 706.005
Accessories	
Aerosol-tight cap*1 for 100 mL rectangular bucket, set of 2	5804 712.005

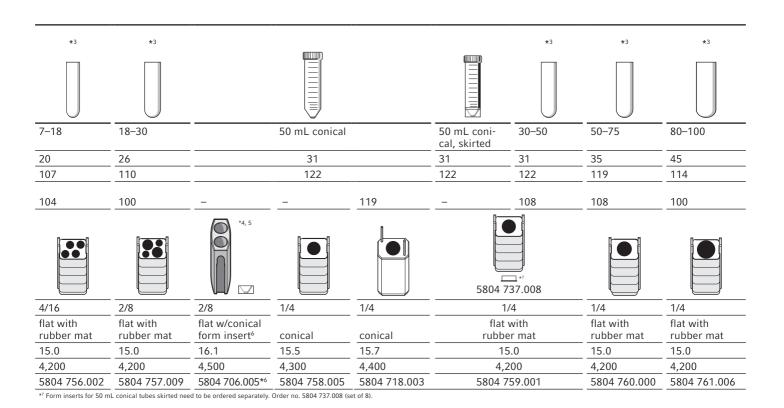

Rotor A-2-DWP-AT and A-2-DWP





Aerosol-tight*1 deepwell plate Rotor A-2-DWP-AT*2

- > For use in Centrifuge 5810/5810 R only
- > Max. speed: 3,486 \times g (4,500 rpm)
- > Rotor can handle up to
 - -2 × 6 MTP
 - -2×2 DWP
 - -2×1 Filtration plate kits
- > Maximum loading height of 67 mm (89 mm without lid)
- > Rotor comes with aerosol-tight caps*1 for safe centrifugation of hazardous samples
- > Rotor and buckets are autoclavable

Deepwell plate Rotor A-2-DWP


- > For use in Centrifuge 5804/5804 R and 5810/5810 R
- > Max. speed: 2,250 \times g (3,700 rpm)
- > Rotor can handle up to
 - $-2 \times 5 MTP$
 - -2×4 Cell culture plates
 - -2×2 DWP
 - -2×1 Filtration plate kits
- > Maximum loading height of 89 mm
- > Rotor and buckets are autoclavable

Description	
Aerosol-tight deepwell plate Rotor A-2-DWP-AT*2	
Rotor A-2-DWP-AT, incl. buckets, aerosol-tight caps and plate carrier	
Deepwell plate Rotor A-2-DWP	
Rotor A-2-DWP, incl. 2 deepwell plate buckets	

^{*1} Aerosol-tightness is tested and certified by Public Health England, Porton Down, UK.

*2 Please note that a software upgrade may be necessary for older centrifuges when using Rotor A-2-DWP-AT.

Please contact your local Eppendorf organization or dealer for further information.

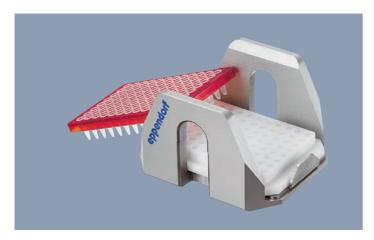




PCR plate adapter (PCR 96)

- > PCR plate adapters are designed to support thin-walled PCR plate wells during centrifugation
- > Available for 96-well and 384-well PCR plates
- > Adapters can be used in all Eppendorf plate rotors

IsoRack adapter

- > IsoRack adapters to support 0.5 mL and 1.5/2 mL IsoRacks during centrifugation
- > Adapters can be used in Eppendorf plate rotors A-2-DWP, A-4-81-MTP/Flex and S-4-104

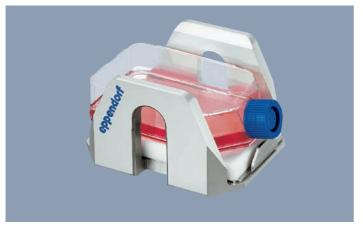



Cell culture flask adapter

- > Adapter to spin down cell cultures directly in a cell culture flask
- > Suitable for 25 cm² flasks
- > Adapters can be used with Rotor A-4-81-MTP/Flex

CombiSlide® adapter

- > Designed for drying slides after hybridization
- > Adapter can hold up to 12 regular glass slides
- > Removable inserts allow centrifugation of glass slides across and parallel to the direction of rotation
- > Adapters can be used in all Eppendorf plate rotors

Order no.
5825 711.009
5825 713.001
5825 719.000
5825 708.008
5825 709.004
5825 706.005

Rotor F-34-6-38




Product features

- > High-speed rotor for molecular applications in tubes from 0.2 mL to 85 mL
- > Max. speed 18,500 \times g (12,000 rpm) in Centrifuge 5810 R or $15,500 \times g$ (11,000 rpm) in 5804/5804R/5810
- > Low rotor weight (3.2 kg) for easy handling
- > 34° borehole angle reduces pellet smears along tube wall
- > Anodized aluminum for high chemical resistance
- > Rotor, lid and adapters are autoclavable

Tube					Δ
Common volumes (mL)	1.5/2	2–5 13 × 75 mm	4–7 13 × 100 mm	4–10	Eppendorf Tubes® 5.0 mL
Max. tube diameter (Ø mm)	11	13	13	16	17
Max. tube height for use lid (mm)	110	90	110	112	104
Adapter options for Rotor F-34-6-38					() priposeth
Capacity per adapter/rotor	4/24	3/18	3/18	2/12	1/6
Adapter bottom shape	round	flat	flat	round	conical
Centrifugation radius (cm)	11.3	10.6	11.4	11.2	10.5
RCF (g-force) at at 12,100 rpm (5810 R) at 11,000 rpm (5804/5804 R/5810)	18,200 15,300	17,070 14,340	18,350 15,420	18,000 15,150	16,824 14,150
Order no. (set of 2)	5804 770.005	5804 738.004	5804 739.000	5804 771.001	5804 777.000
*1 Max. speed in Centrifuge 5810 R, lower speed applies to					

^{*2} Nominal tube volume as published by the tube manufacturer.

Description	Order no.
Rotor F-34-6-38 for 6 × 85 mL tubes incl. rotor lid	5804 727.002

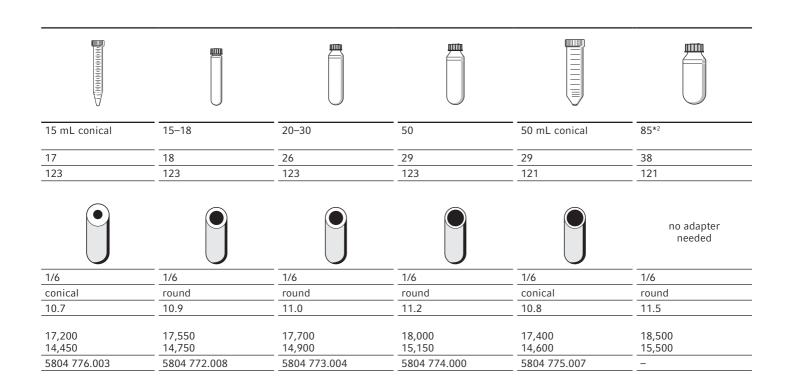

Rotor FA-45-6-30



Product features

- > High-speed rotor for applications in tubes from 5–50 mL
- > Max. speed 20,130 \times g (12,100 rpm) in Centrifuge 5810 R or 16,639 $\times g$ (11,000 rpm) in 5804/5804R/5810
- > Eppendorf QuickLock rotor lid for fast lid locking
- > Aerosol-tight*1 for safe centrifugation of hazardous samples
- > Low rotor weight (3.2 kg) for easy lifting
- > Anodized aluminum for high chemical resistance
- > 45° borehole angle minimizes pellets smear
- > Rotor, lid and adapters are autoclavable

•	Description	Order no.
	Rotor FA-45-6-30*3, incl. aerosol-tight Eppendorf QuickLock® lid	5820 715.006


Tube					A
Common volumes (mL)	2.6–5 13 × 75 mm	4–8 13 × 100 mm	5.5–10 16 × 75 mm	7.5–12 16 × 100 mm	Eppendorf Tubes® 5.0 mL
Max. tube diameter (Ø mm)	13	13	16	16	17
Max. tube height for use with aerosol-tight lid*1 (mm)	119	119	119	119	121
Adapter options for Rotor FA-45-6-30	O propuseddo			Juspuneddie	
Capacity per adapter/rotor	1/6	1/6	1/6	1/6	1/6
Adapter bottom shape	round	round	round	round	conical
Centrifugation radius (cm)	12.0	12.0	12.0	12.0	12.1
RCF (g-force) at at 12,100 rpm (5810 R) at 11,000 rpm (5804/5804 R/5810)	19,246 16,233	19,246 16,233	19,246 16,233	19,246 16,233	19,806 16,369
Order no. (set of 2)	5820 726.008	5820 725.001	5820 728.000	5820 727.004	5820 730.005

^{*1} Aerosol-tightness tested by Public Health England, Porton Down, UK.
*2 Max. speed in Centrifuge 5810 R, lower speed applies to Centrifuge 5804/5804R/5810.
*2 Please note that a software upgrade may be necessary for older centrifuges when using Rotor FA-45-6-30.
Please contact your local Eppendorf organization or dealer for further information.

The New **Eppendorf Conical Tubes**

Whether centrifugation, mixing, or other applications the Eppendorf Conical Tubes 15 mL and 50 mL will perfectly fit into your workflows. They represent the next step in Eppendorf's comprehensive system solutions in liquid, sample and cell handling.

More information: www.eppendorf.com/conicals

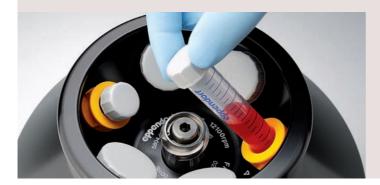



7–17	15 mL conical	15–18	20–30	35	50 mL conical
17.5	17	18	26	29	31
122	125	125	125	125	121
propuedibe		paguedde	Drapuneddio		no adapter needed
1/6	1/6	1/6	1/6	1/6	1/6
1/6 round	conical	1/6 round	round	conical	conical
1/6 round		1/6			
1/6 round 12.0 19,246 16,233	conical	1/6 round	round	conical	conical

Rotor F-35-48-17 and FA-45-20-17





Rotor F-35-48-17*1

- > High-capacity fixed-angle rotor for 15 mL conical and glass tubes
- > Max. speed: 5,005 \times g (5,500 rpm)
- > Form inserts for 15 mL conical tubes

Rotor FA-45-20-17*1

- > 20 places for Eppendorf Tubes® 5.0 mL
- > Max. speed: 20,913 \times g (13,100 rpm)
- > Eppendorf QuickLock rotor lid for fast lid locking
- > Rotor, lid and adapters are autoclavable
- > Anodized aluminum for high chemical resistance

Description	Order no.
Fixed-angle Rotor F-35-48-17	
Includes 48 steel sleeves for 15 mL tubes, with 40 conical inserts and 48 rubber mats	5820 772.000
Includes 24 steel sleeves for 15 mL tubes, with 24 conical inserts and 24 rubber mats	5820 771.003
Fixed-angle rotor FA-45-20-17	
Rotor FA-45-20-17, 20 × 5.0 mL, incl. aerosol-tight Eppendorf QuickLock® rotor lid	5820 765.003
Adapter for 1.5/2.0 mL tubes, set of 4	5820 768.002
Adapter for cryogenic vials, set of 4	5820 769.009
Adapter for HPLC tubes, set of 4	5820 770.007

^{*}¹ Please note that a software upgrade will be necessary for older centrifuges when using Rotor F-35-48-17 or FA-45-20-17. Please contact your local Eppendorf organization or dealer for further information.





Rotors F-45-30-11 and FA-45-30-11

- > 30 places for 1.5/2.0 mL tubes
- > Max. speed: 20,800 \times g (14,000 rpm)
- > Rotor FA-45-30-11 comes with aerosol-tight*2 lid for safe centrifugation of hazardous samples
- > Rotors and lids are autoclavable

Rotor FA-45-48-11*3

- > 48 places for 1.5/2.0 mL tubes
- > Max. speed: 19,083 \times g (max. 13,000 rpm)
- > Eppendorf QuickLock rotor lid for fast lid locking
- > Aerosol-tight*2 for safe centrifugation of hazardous samples
- > Anodized aluminum for high chemical resistance
- > Rotor and lid are autoclavable

Drum Rotor T-60-11

- > High capacity for $60 \times 1.5/2.0$ mL tubes
- > Max. speed: 16,400 \times g (14,000 rpm)
- > Horizontal centrifugation for pellet formation at tube bottom—ideal for low sample quantities

Rotor F-45-48-PCR

- > Accommodates up to 6 \times 5/8-tube PCR strips or 48×0.2 mL PCR tubes
- > Safe and easy handling without additional lid
- > Max. speed: 15,350 $\times q$ (12,000 rpm)



Description	Order no.
Rotor FA-45-48-11, incl. aerosol-tight Eppendorf QuickLock® rotor lid	5820 760.001
Rotor FA-45-30-11, 30 × 1.5/2.0 mL, with aluminum lid, aerosol-tight*1	5804 726.006
Rotor F-45-30-11, 30 × 1.5/2.0 mL, with aluminum lid	5804 715.004
Adapter for 0.2 mL PCR tubes (for 1.5/2 mL rotors), set of 6	5425 715.005
Adapter for 0.4 mL tubes (for 1.5/2 mL rotors), set of 6	5425 717.008
Adapter for 0.5 mL tubes (for 1.5/2 mL rotors), set of 6	5425 716.001
Adapter for BD Microtainer® (for 1.5/2 mL rotors), set of 6	5425 716.001
Drum rotor T-60-11, with lid, without adapters	5804 730.003
Adapter for 10 × 1.5/2 mL tubes (Ø 11 mm), set of 6	5804 731.000
Adapter for 20 × 0.4 mL tubes (Ø 6 mm), set of 6	5804 732.006
Rotor F-45-48-PCR, for 6 × 8-tube PCR strips	5804 735.005


 ^{*2} Aerosol-tightness is tested and certified by Public Health England, Porton Down, UK.
 *3 Please note that a software upgrade will be necessary for older centrifuges when using Rotor FA-45-48-11.
 Please contact your local Eppendorf organization or dealer for further information.

Rolling Cabinet




If space is a limited resource in your lab, Eppendorf's rolling cabinet can give you the freedom and flexibility to use your Centrifuge 5804/5804 R or 5810/5810 R wherever you need it. Use the rolling cabinet if there is not enough space on the regular lab bench or to transport and use the centrifuge in other labs.

The included drawer provides extra storage space for your rotors and adapters.

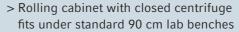



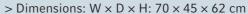



Watch video!

Ordering information

Description	Order no.
Rolling cabinet, for Centrifuge 5804/5804 R and 5810/5810 R	5821 020.010

Centrifuge 5804/5804 R and 5810/5810 R

Technical specifications

Description	Centrifuge 5804	Centrifuge 5804 R	Centrifuge 5810	Centrifuge 5810 R			
Max. rcf							
> Fixed-angle rotor	20,913 × g	20,913 × <i>g</i>	20,913 × g	20,913 × g			
> Swing-bucket rotor	4,500 × <i>g</i>	4,500 × <i>g</i>	4,500 × <i>g</i>	4,500 × g			
> Plate rotor	2,250 × g	2,250 × <i>g</i>	3,486 × <i>g</i>	3,486 × g			
Speed		200–14,000 rpm (rpm (in 10 rpm increments)				
Max. capacity	4 × 250 mL	4 × 250 mL	4 × 750 mL	4 × 750 mL			
Rotor options	12	12	18	18			
Accel./Braking ramps		10/10					
Programs	35 user programs						
Timer	1 min. to 99 min, with continuous mode, short spin						
Temperature control range	N/A	N/A -9 °C to 40 °C N/A		−9 °C to 40 °C			
Noise level							
> Rotor S-4-104 (4 × 750 mL)			<70 dB(A)	<56 dB(A)			
> Rotor S-4-72 (4 × 250 mL)	<62 dB(A)	<58 dB(A)					
> Rotor FA-45-6-30 (6 × 50 mL)	<55 dB(A)	<54 dB(A)	<56 dB(A)	<55 dB(A)			
Power supply	230 V/50–60 Hz			230 V/50-60 Hz			
Power requirement	Max. 900 W	Max. 1,650 W	Max. 900 W	Max. 1,650 W			
Dimensions (W × D × H)	47 × 55 × 34 cm	64 × 55 × 34 cm	54 × 61 × 35 cm	70 × 61 × 35 cm			
Access height	28 cm	28 cm	28 cm	28 cm			
Height with open lid	74 cm	74 cm	80 cm	80 cm			
Weight without rotor	 55 kg	80 kg	68 kg	99 kg			

Centrifuge Packages



Ordering	information
Oraciniq	minormation

Cell culture centrifuge packages*	Max. capacity of conical tubes per rotor	Order no.
Centrifuge 5804, with Rotor A-4-44 incl. adapters for 15 mL/50 mL conical tubes	16 × 15 mL / 4 × 50 mL	5804 000.323
Centrifuge 5804 R, refrigerated, with Rotor A-4-44 incl. adapters for 15 mL/50 mL conical tubes	16 × 15 mL / 4 × 50 mL	5805 000.327
Centrifuge 5804, with Rotor S-4-72 incl. adapters for 15 mL/50 mL conical tubes	32 × 15 mL / 16 × 50 mL	5804 000.528
Centrifuge 5804 R, refrigerated, with Rotor S-4-72 incl. adapters for 15 mL/50 mL conical tubes	32 × 15 mL / 16 × 50 mL	5805 000.629
Centrifuge 5810, with Rotor A-4-62 incl. adapters for 15 mL/50 mL conical tubes	36 × 15 mL / 16 × 50 mL	5810 000.327
Centrifuge 5810 R, refrigerated, with Rotor A-4-62 incl. adapters for 15 mL/50 mL conical tubes	36 × 15 mL / 16 × 50 mL	5811 000.320
Centrifuge 5810, with Rotor A-4-81 incl. adapters for 15 mL/50 mL conical tubes	48 × 15 mL / 20 × 50 mL	5810 000.424
Centrifuge 5810 R, refrigerated, with Rotor A-4-81 incl. adapters for 15 mL/50 mL conical tubes	48 × 15 mL / 20 × 50 mL	5811 000.428
Centrifuge 5810, with Rotor S-4-104 incl. adapters for 15 mL/50 mL conical tubes	56 × 15 mL / 28 × 50 mL	5810 000.521
Centrifuge 5810 R, refrigerated, with Rotor S-4-104 incl. adapters for 15 mL/50 mL conical tubes	56 × 15 mL / 28 × 50 mL	5811 000.622
Centrifuge packages for round-bottom tubes	Max. capacity of round- bottom tubes per rotor	Order no.
Centrifuge 5810, with Rotor S-4-104 incl. adapters for round-bottom tubes with diameter 13/16 mm NEW	92 × 13 mm / 80 × 16 mm	5810 000.629
Centrifuge 5810 R, refrigerated, with Rotor S-4-104 incl. adapters for round-bottom tubes with diameter 13/16 mm NEW	92 × 13 mm / 80 × 16 mm	5811 000.827
ADDRIVED ON I		• ———

^{*230} V/50-60 Hz Technical specifications subject to change. Errors and omissions excepted.

eppendorf

Eppendorf Microcentrifuges



















									-
	MiniSpin®	MiniSpin plus	5418	5418 R	5424	5424 R	5427 R	5430	5430 R
Model	Entry-level	Entry-level		apacity for ar biology	The laborat	tory standard	High throughput		rifuge with e capabilities
Capacity	12 × 1.5/2 mL	12 × 1.5/2 mL	18 × 1.	.5/2 mL	24 × 1	.5/2 mL	48 × 1.5/2 mL		5/2 mL nL, 2 × MTP)
Max. speed (RCF)	12,100 × g	14,100 × g	16,87	73 × g	20,238 × g	21,130 × g	25,000 × g	30,13	30 × g
Max. rpm	13,400	14,500	14,	.000	14,680	15,000	16,220	17,	500
Rotor options	2	2		1		4	9	1	2
Temperature settings				0 °C to 40 °C		-10 °C to 40 °C	-10 °C to 40 °C		-11 °C to 40 °C
Order no.	5452 000.018	5453 000.011	5418 000.017	5401 000.013	5424 000.614	5404 000.618	5409 000.012	5427 000.615	5428 000.619

Eppendorf Multipurpose Centrifuges















	5702	5702 R	5702 RH	5804	5804 R	5810	5810 R
Model	medium capacity	fuge optimized for y needs in clinical cell culture labs	Active heating and cooling for high temperature accuracy to maximize cell viability (e.g. cell transfection)	Compact high speed centrifuge for medium capacity needs. With swing-bucket rotor and deepwell plate capacity		The workhorse for mid to high-throughput laboratories: Highly versatile, high speed and high plate capacity	
Capacity	4 × 100 mL			4 × 250 mL		4 × 750 mL	
- Conical tubes	20 × 15 mL/4 × 50 ml		ηL	40 × 15 mL/16 × 50 mL		56 × 15 mL/28 × 50 mL	
- Blood collection tubes	36 × 13 mm/30 × 16 mm		mm	56 × 13 mm/52 × 16 mm		100 × 13 mm/80 × 16 mm	
– Plates				10 × MTP/2 × DWP		16 × MTP/4 × DWP	
Max. speed (RCF)	3,000 × g			20,913 × g		20,913 × g	
Max. rpm	4,400			14,000		14,000	
Rotor options	6			12		18	
Temperature settings		−9 °C to 40 °C	−9 °C to 42 °C		−9 °C to 40 °C		−9 °C to 40 °C
Order no.	5702 000.019	5703 000.012	5704 000.016	5804 000.013	5805 000.017	5810 000.017	5811 000.010

Technical specifications subject to change. Errors and omissions excepted.

Your local distributor: www.eppendorf.com/contact

Eppendorf AG 22331 · Hamburg · Germany E-mail: eppendorf@eppendorf.com · www.eppendorf.com

www.eppendorf.com/centrifugation

BD Vacutainer® and BD Falcon® are registered trademarks of Becton, Dickinson and Company, Microtainer® is a registered trademark of Becton Dickinson, New Jersey, USA. Centriprep® is a registered trademark of Millipore Corporation. Corning® is a registered trademark of Corning, Inc. DURAN® is a registered trademark of DURAN GROUP GmbH, Germany. Eppendorf®, the Eppendorf logo, Eppendorf DuckLock®, Eppendorf Tubes®, CombiSlide®, epGreen®, epServices® logo and MiniSpin® are registered trademarks of Eppendorf AG, Germany. U.S. Design Patents are listed on www.eppendorf.com/jp - All rights reserved, including graphics and images. Copyright © 2014 by Eppendorf AG. Order No. A58X X12 020/GB8/8T/1214/CCHH/KGR · Carbon neutrally printed in Germany.