## IKA® RET control-visc

Betriebsanleitung	DE	4
Ursprungssprache		
Operating instructions	EN	20
Mode d'emploi	FR	36
Руководство пользователя	RU	52
使用说明	ZH	68
取扱説明書	JA	84
사용 설명서	KO	100

Indicaciones de seguridad	ES	116	Biztonsági utasítások	HU	126
Veiligheidsinstructies	NL	117	Varnostna opozorila	SL	127
Norme di sicurezza	IT	118	Bezpečnostné pokyny	SK	128
Säkerhetsanvisningar	SV	119	Ohutusjuhised	ET	129
Sikkerhedshenvisninger	DA	120	Drošības norādes	LV	130
Sikkerhedsanvisninger	NO	121	Nurodymai dėl saugumo	LT	131
Turvallisuusohjeet	FI	122	Инструкции за безопасност	BG	132
Instruções de serviço	PT	123	Indicații de siguranță	RO	133
Wskazówki bezpieczeństwa	PL	124	Υποδείξεις ασφάλειας	EL	134
Bezpečnostní pokyny	CS	125			

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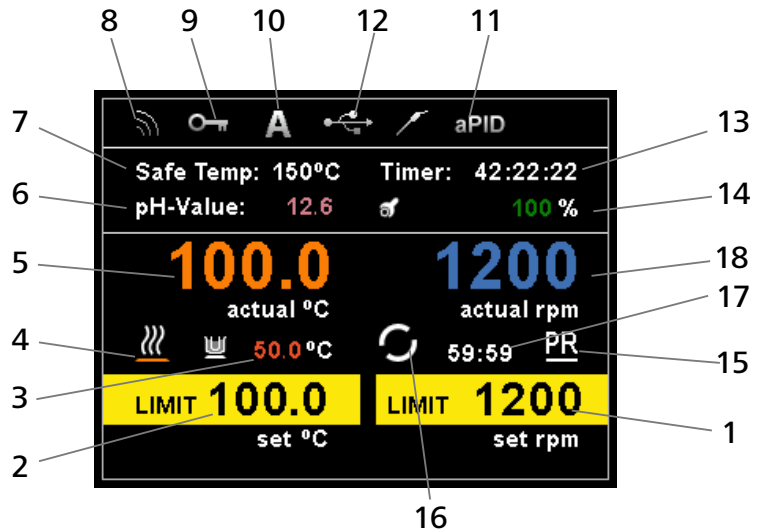
- 1 Set value speed
- 2 Set value temperature
- 3 Temperature limitation of carrier liquid  
(only in connection with double temp-sensors)
- 4 Heater activated
- 5 Actual temperature
- 6 Measured pH-Value (only in connection with pH-probe)
- 7 Actual safety circuit temperature
- 8 Wireless connection to remote control
- 9 All operation elements locked
- 10 Operation mode (A/B/C)
- 11 External temperature probe connected  
and the selected control mode
- 12 USB connection with PC
- 13 Timer / Counter
- 14 Viscosity trend value
- 15 Connection with PC established
- 16 Motor activated
- 17 Intermittent status
- 18 Actual speed value

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## Contents

	Page		
Controls and display	2	Menu(overview)	25
Declaration of Conformity	20	Connections and interfaces	30
Warranty	20	Accessories	32
Explication of warning symbols	20	Troubleshooting	33
Safety instructions	21	Maintenance and cleaning	34
Unpacking	22	Technical data	34
Correct use	22		
Assembling the support rod	22		
Setting the safety temperature limit	23		
Operation	24		

## Declaration of Conformity

We declare under our sole responsibility that this product is in compliance with the regulations 2014/35/EU, 2014/30/EU and 2011/65/EU and conforms to the standards or normative documents: EN 61010-1, EN 61010-2-010, EN 61010-2-051, EN 61326-1, EN 60529 and EN ISO 12100.

## Warranty

According to **IKA**®'s Terms and Conditions of sale and delivery, this product is covered by a warranty for a period of 24 months. In case of making a warranty claim, please contact your local dealer or, if you wish, you can send the device directly to our factory. Please include the sales invoice and state the reasons for your guarantee claim. In this case, you are responsible for shipping and handling costs.

The warranty does not cover wearing parts, nor to defects that are the result of improper use, insufficient care and maintenance or failure to follow the instructions in this operating manual.

## Explication of warning symbols

General hazard

**DANGER**

This symbol indicates information which is **essential for your health**. Failure to observe this information may be detrimental to health or may result in injuries.

**WARNING**

This symbol indicates information which is important for **ensuring that the device functions without any technical problems**.

Failure to observe this information could damage the device.

**CAUTION**

This symbol indicates information **which is important for proper use of the device and/or ensuring that the device functions correctly**. Failure to observe this information can lead to inaccurate results.

**ATTENTION**

ATTENTION - risk of damage due to magnetism.

**DANGER**

DANGER - note on hazards arising from a hot surface.

## Safety instructions

- **Please read the instruction manual in full before use and follow the safety instructions.**

- Keep the instruction manual in a place where it can be accessed easily.

### General instructions

- Ensure that only trained staff use the device.
- Be sure to comply with all safety instructions, directives and all matters of health, safety and accident prevention in the workplace.
- The plug must be earthed (grounded contact).

**ATTENTION**

### Caution - magnetism!

Beware of possible effects from the magnetic field (pacemakers, data media, etc.).

**DANGER**

### Burns hazard!

Exercise caution when touching housing components and the heating plate. The

heating plate can reach temperatures of up to 320 °C. Watch for residual heat after appliance has been switched off.

**The device should only be transported once it has cooled!**

### Device configuration

**DANGER**

Do **not** operate the device in explosive atmospheres, in the presence of hazardous materials or under water.

- Install the device on a flat, stable, clean, non-slip, dry and fire-resistant surface.
- The device feet must be clean and undamaged.
- The power supply cable and cables to the external sensors must not be allowed to come into contact with the heated mounting plate.
- Prior to each use, always check the device and accessories for damage. Never use damaged parts.

### Permissible media / Contamination / Side reactions

**WARNING**

**Caution!** Only media whose flashpoint lies above the set safety temperature limit (50 °C - 380 °C) may be processed or

heated with this device.

The set safety temperature limit must be at least 25 °C below the flashpoint of the medium in use.

**WARNING**

Pay attention to the risks associated with:

- flammable materials
- flammable media with low steam pressure
- glass breakage
- incorrect vessel size
- overfilling of medium
- unstable vessel.
- Only process pathogenic materials in closed vessels under a suitable extractor hood. Please contact **IKA®** if you have any questions.

**DANGER**

Only work with media for which the energy input during processing is harmless. This also needs to take into account other

sources of energy such as, for example, light irradiation.

- At high speeds the mounting plate is liable to warming, even if the device is not in heating mode.
- Bear in mind the possibility of contamination or unwanted chemical reactions.
- There is the possibility that friction from rotating accessory parts can result in contamination of the medium being processed.
- Bear in mind the following when using PTFE-coated magnetic rods: *Chemical reactions of PTFE occur in contact with molten or solute alkali metals and alkaline earth metals as well as with fine powders of metals in groups 2 and 3 of the periodic table*

at temperatures above 300 - 400 °C. Only elementary fluorine, chlorine trifluoride and alkali metals attack it; halogenated hydrocarbons have a reversible swelling effect.

(Source: Römpf "Chemie-Lexikon" and "Ullmann" Vol. 19)

- Only glass-coated magnetic rods should be used in conjunction with solute alkali metals or alkaline earth metals or at temperatures above 250 °C.

### Performing trials

**WARNING**

Always wear personal protective equipment in accordance with the hazard class of the media being worked with.

Otherwise there is the risk of the following:

- spraying and evaporation of liquids
- ejection of parts
- release of toxic or inflammable gases.
- If any of the following happens, reduce the speed:
  - medium sprays out of vessel due to excessive speed
  - process is not smooth
  - vessel moves around on the mounting plate.

### Accessories

- Safe operation can only be ensured when working with accessories as described in the "**Accessories**" section.
- Only put together accessories with the power supply cable disconnected.
- Refer to the operating instructions for the accessories.
- Ensure that the external temperature sensor is inserted into the medium to a depth of at least 20 mm when connected.
- Accessories must be connected securely to the device and must not come loose on their own. The centre of gravity of the device must be located within the mounting plate.

### Power supply / switching off the device

- The specified settings on the rating plate must coincide with the actual power supply.
- The plug for the power supply connection must be easy to access and remove.
- The device is only disconnected from the power supply when the power or device plug is removed.

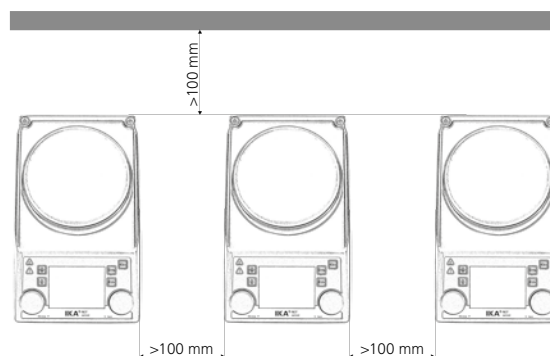
**WARNING**

After an interruption to the power supply, the device starts up again in Mode B.

### Protection of the device

- The device must only be opened by trained, skilled personnel.
- Do not cover the device, even partially, with elements such as metallic plates or sheets otherwise it may overheat.
- Avoid knocks or impact to the device or accessories.
- Ensure the mounting plate remains clean.
- Observe minimum distances between devices. Between device and wall should be 100 mm (min), above the assembly should be 800 mm (min).

## Unpacking

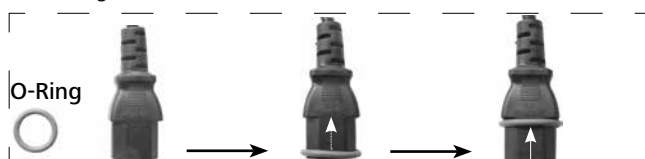
### • Unpacking

- Unpack the device carefully
- Any damage should be notified immediately to the shipping agent (post office, railway network or transport company)

### • Scope of delivery

- RET control-visc heated magnetic stirrer
- Power supply cable
- Operating instructions
- Screwdriver (safety circuit)
- USB cable

- Protective cover: **H 104**
- Magnetic stirring bar: **IKAFLON 30** and **40** mm
- Temperature sensor: **PT 100.70**
- O Ring

Please slide the O-Ring on the power cord plug to avoid ingress of water into the connector.

## Correct use

### • Use

- for mixing and/or heating liquids.

### • Area of use (only indoors)

- laboratories
- pharmacies
- schools
- universities

The device is suitable for use in all areas apart from domestic areas and areas directly connected to a low-voltage supply which also serves domestic areas.

The safety of the user cannot be guaranteed if the device is used in conjunction with accessories which are not supplied or recommended by the manufacturer, if the device is used contrary to the intended purpose against the manufacturer's instructions or if the device itself or PCB has been subjected to modifications by third parties.

## Assembling the support rod

- Remove the threaded plug (**P**)
- Remove the protective cap from the support rod
- Place the washer between the housing and the nut
- Screw in the support rod manually until it cannot be tightened further
- Tighten the nut using an SW17 open-end spanner
- Assemble the accessories using a cross sleeve

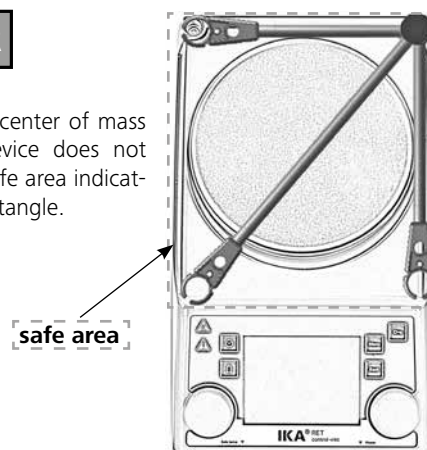
### Note:

When using bath attachments with a diameter over 180 mm, please use support rod in conjunction with an extension. (See "Accessories")

### Risk of tipping!

Make sure that the center of mass of the attached device does not protrude over the safe area indicated with a dotted rectangle.

## Setting the safety temperature limit

The maximum achievable heating plate temperature is restricted by an adjustable safety temperature limit. Once this limit has been reached, the device stops heating.

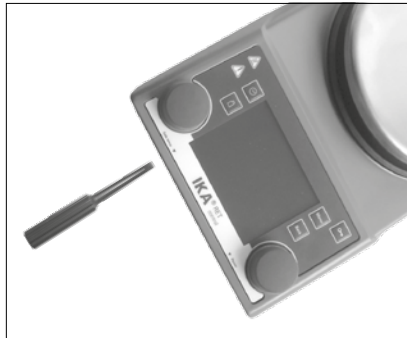
**WARNING**

**The safe temperature limit must always be set at least 25 °C lower than the flash point of the media to be processed!**

The adjustable maximum heating plate temperature is at least 10 °C under the set safety temperature limit.

Setting range: [50 °C] to [max. SET temperature + 50 °C]

Factory setting: [max. SET temperature + 50 °C]

### Setting the safety temperature limit

After switching on the device, the safety temperature limit (7) can be adjusted using a screwdriver.

Do not turn the setting screw beyond the clockwise or anticlockwise stop. This will cause irreparable damage to the potentiometer.

- Set the main switch (A) to the ON position
- Using the screwdriver supplied, turn the "Safe Temp" setting screw (B) to the clockwise stop.
- Use the temperature rotary knob (C) to set the target temperature to the desired "Safe Temperature" and wait until the temperature is achieved.
- Turn the "Safe Temperature" setting screw (B) slowly anticlockwise until the heating function switches off and the display shows [Er 25].
- The "Safe Temp" value (7) is displayed on the screen.

**WARNING**

Setting the safety temperature limit as to be done as described.

The value "Safe Temperature" shown on the display only serves for visualization.

### Functional test: safety circuit switch-off

- Heat the device to a temperature of over 50 °C.
- Set the safety temperature limit as far to the left as possible (50 °C) and set the mains switch (A) to the ON position.
- The display will show: [Er 25]

## Operation

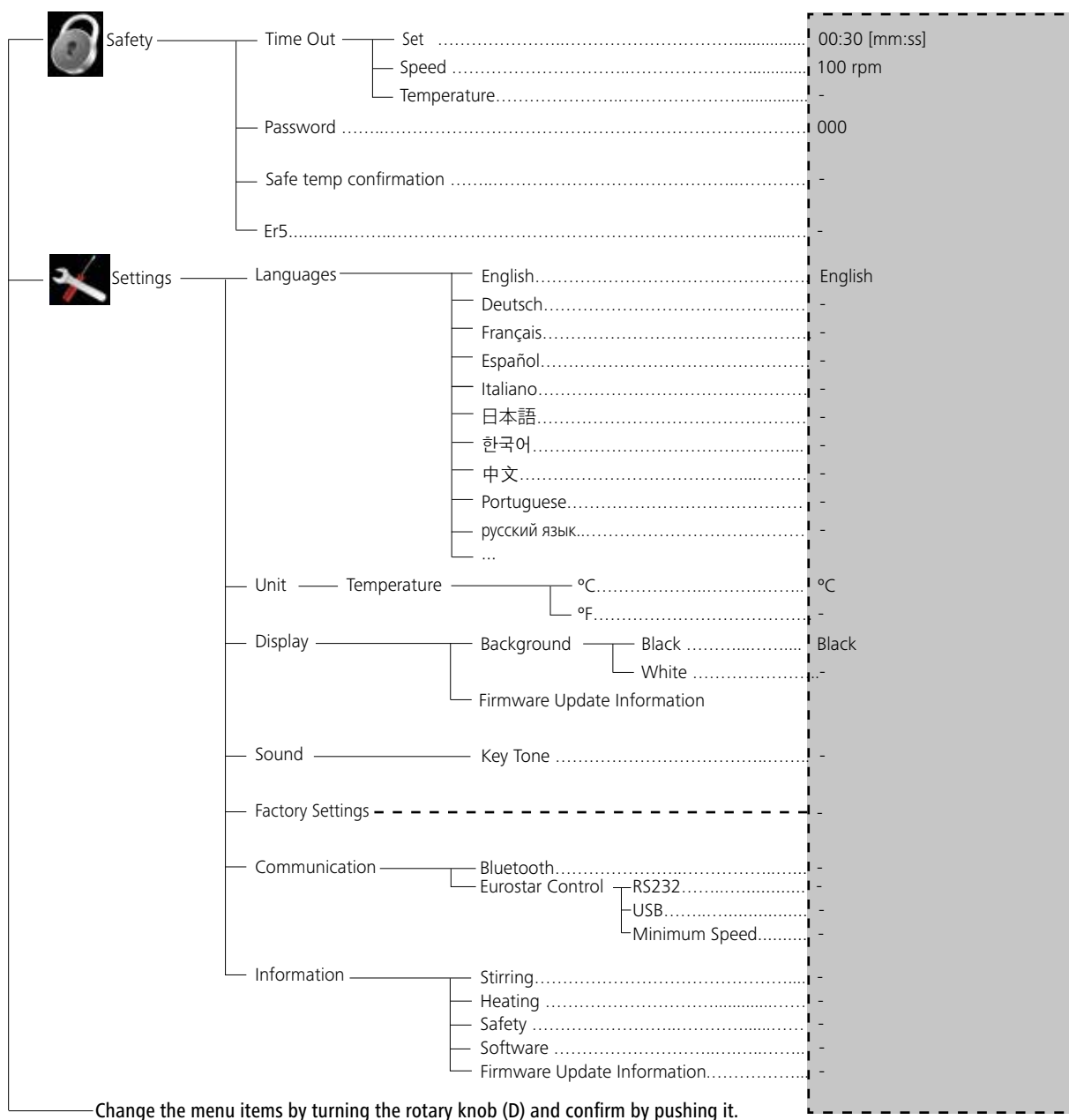
Remove the protective film from the mounting plate before use! The "Back" menu key is used to change between menu items.

<b>Start-up</b>	<ul style="list-style-type: none"> <li>☞ Set the main switch (A) to the OFF position</li> <li>☞ Insert the power supply cable into the mains socket (K)</li> <li>☞ Set the main switch (A) to the ON position</li> <li>⇒ Standard mode A is selected (see "Menu (details)")</li> </ul>
<b>Stirring</b>	<ul style="list-style-type: none"> <li>☞ Set the speed using the rotary knob (D)</li> <li>⇒ The set value is shown on the display (J)</li> <li>☞ Press the rotary knob (D) to start the stirring function</li> </ul> <p>① ⇒ The symbol (16) appears when the motor is active</p>
<b>Heating</b>	<ul style="list-style-type: none"> <li>☞ Set the safety temperature limit (see "Menu (details)")</li> <li>☞ Set the target temperature using the rotary knob (C)</li> <li>⇒ The set value is shown on the display (J)</li> <li>☞ Set the temperature control mode (see "Menu (details)")</li> <li>☞ Press the rotary knob (C) to start the heating function</li> </ul> <p style="text-align: center;">The target and current temperature are shown permanently on the display (J):</p> <p>① ⇒ When the heating function is switched on, the heating symbol (4) is displayed.</p> <div style="border: 1px solid black; padding: 2px; width: fit-content; margin: 5px auto;">The maximum adjustable heating plate temperature setting is 340 °C</div> <p>① ⇒ Whenever the mounting plate is at a temperature above 50 °C, the display (J) shows  and the current temperature is shown, even if the device is switched off.</p>
<b>Connecting an external thermometer</b> <i>(direct temperature control in the medium)</i>	<ul style="list-style-type: none"> <li>☞ Set the main switch (A) to the OFF position</li> <li>☞ Connect the safety contact thermometer compliant with DIN 12878 Class 2 or temperature sensor PT 100 (individual sensor) or PT 1000 (double sensor) to socket (L)</li> <li>☞ Set the main switch (A) to the ON position</li> </ul> <p>①                      Temperature sensor PT 100    ⇒ The current temperature (5) for the temperature sensor shown on the display (J) corresponds to the temperature of the medium. The "External temperature probe" symbol (11) appears.</p> <p>①                      Temperature sensor PT 1000    ⇒ The current temperature (5) for the temperature sensor shown on the display (J) corresponds to the temperature of the medium. The "External temperature probe" symbol (11) appears. The temperature of the carrier fluid (3) is also displayed.</p>
<b>Weighing</b>	<ul style="list-style-type: none"> <li>☞ Press the (G) key</li> <li>☞ Wait until the system is stable (once the system is stable, a "g" appears after the digit 0).</li> <li>☞ Place the weight in the centre of the heating plate and wait until the system is stable again.</li> </ul> <p>① ⇒ The maximum display range is up to 2000 g</p> <p>⇒ The weighing range is max. 5000 g</p> <p>⇒ It is only possible to use the weighing mode when the heating or stirring functions are not active.</p> <p>① ⇒ The weighing module can be calibrated.</p>
<b>pH measurement</b>	<ul style="list-style-type: none"> <li>⇒ See section entitled "Menu (details)"</li> </ul>
<b>Timer</b>	<ul style="list-style-type: none"> <li>☞ Press the (H) key</li> <li>☞ Set the time in hours, minutes and seconds by turning the knob (D) and confirm by pressing it; the timer function starts to count down the remaining time once the heating or stirring function has started (see "Menu (details)").</li> </ul> <p>①    Once the timer reaches 0, the heating function is switched off, even if the motor has only just been started. If the heating is activated during the timer sequence, only the heating is switched off, but the motor continues to run (delay in boiling).</p>

# Menu (overview)

## Menu structure

		Factory settings			
Menu	 Stirring	Torque trend measurement.....	-		
		Stirring bar decoupling.....	-		
		Intermittent Mode	Run/Stop .....	-	
			Interval	Run Time.....	00:00 [mm:ss]
				Stop Time.....	00:00 [mm:ss]
	Speed limit.....	1700 rpm			
	 Heating	Control method	Fast PID (fPID).....	fPID	
			Accurate PID (aPID).....	-	
			Two - point (2pt).....	-	
		Limits	Medium limit.....	-	
			Carrier limit.....	-	
			Limit information.....	-	
	Temperature sensor	Calibration	2-point Calibration .....	-	
			3-point Calibration .....	-	
		Reset Calibration.....	-		
	 Weighing	Measurement .....	-		
		Calibration	2-point Calibration.....	200...5000 g	
			3-point Calibration.....	200...5000 g	
	 Timer	Set.....	00:00:00 [hh:mm:ss]		
		Display.....	-		
 pH probe	Calibration	2-point Calibration.....	-		
		3-point Calibration.....	-		
	Display.....	-			
 Mode	A.....	A			
	B.....	-			
	C.....	-			
 Display	Torque trend measurement.....	-			
	pH value .....	-			
	Timer .....	-			
	Graph .....	-			
 Graph	Axis Assignment	Temperature.....	-		
		Speed.....	-		
		Torque.....	-		
	Axis Scaling	Automatic.....	-		
		Manual.....	-		
 Programs	Program 1.....	-			
	Program 2.....	-			
	Program 3.....	-			
	Program 4.....	-			
	Program 5.....	-			

**Menu (details)**

**Stirring**

Torque trend measurement

The torque trend measurement is used to deduce the change in viscosity of the reaction medium.

The devices are **not** designed to measure absolute viscosity. They only measure and display the relative change in the viscosity of the medium from a starting point specified by the user.

Once the speed of the motor and magnetic rod in the medium have stabilised to the target speed, the viscosity measurement is started at 100%.

Consequently, an increase in the measured torque trend is indicated by a value > 100% and a reduction in values < 100%.

The value can always be reset to 100% by pressing the "Back" (F) membrane key.

**Note:**

**Torque trend measurement only works for a constant set speed for the duration of the measurement.**

**As a result, intermittent mode cannot be used in conjunction with torque trend measurement.**

The current control variable is saved as the reference 100% ΔP and shown on the digital display. The change in the viscosity is then shown in %. Depending on whether the viscosity increases or decreases, the percentage rises or falls above or below 100% respectively.

Stirring bar decoupling

If the device detects that a stirring bar has broken away, the stirring function is interrupted temporarily in order to pick up the bar. The system then returns to the previously set speed. When the system detects a bar decoupling, this is indicated with a special symbol (✘ instead of symbol 16). If another stirring bar decoupling occurs within the next 3 minutes (as long as the target speed has not been modified), once the bar is recovered the system only goes up to a speed of 100 rpm. In the case of 4 consecutive stirring bar decouplings -with less than 3 minutes' normal speed between them - the error message ER 41 is displayed. The speed is reduced to 200 rpm and the heating is switched off.

In such cases, the testing conditions will need to be reconsidered.

**Note:**

The user is responsible for ensuring the suitability/function of the stirring bar decoupling monitoring for the current conditions: stirring bar, speed, vessel, medium, etc.

**Stirring bar decoupling monitoring does not work if the stirring bar already breaks away before the system has reached the set target speed!**

The higher the torque to be transferred, the more safely the stirring bar decoupling monitoring works.

In a glass beaker with water, the stirring bar decoupling for a bar from 40 mm (d=8 mm) upwards can be seen from 300 - 1200 rpm. Longer stirring bars, larger bar surface areas, e.g. Trika, or higher speeds increase the torque to be transferred.

**Heating**

Regulation of the temperature of the medium using PT 100 temperature sensor

When using the PT 100, the selected temperature control mode is shown on the digital display by the  symbol.

The set value temperature (2) and the actual temperature (5) shown on the display refer to the temperature of the medium. The heating plate temperature is limited by the actual safety circuit temperature (7). The medium temperature is controlled by means of a microcontroller. The medium temperature is measured by the PT 100 temperature sensor and is heated up as quickly as possible without overshooting to the set temperature. The microcontroller is capable of adjusting automatically to the different heating capacities of various vessels and their contents. As a result, the heat is distributed in an optimum fashion with no temperature drift or ripples.

Regulation of the temperature of the medium using PT 1000 double temperature sensor

Operation with the PT 1000 double sensor is indicated on the digital display by the  symbol. One temperature sensor is located in the heat carrier and the other is located in the reaction medium. The set value temperature (2) refers to the temperature of the medium. The heat carrier temperature is shown on the display as an additional current temperature with the symbol .

The actual temperature of the reaction medium is shown as the actual temperature (5). The heating plate temperature (and therefore the heat carrier temperature) are limited by the actual safety circuit temperature (7). The reaction medium temperature is limited by the set temperature limits (carrier/reaction temperature limit).

Intermittent mode

Intermittent mode consists of regular, cyclical interruption of the stirring function. The run time and interval can both be set. When switching on the device, the time for each is displayed if intermittent mode is activated.

Speed limit

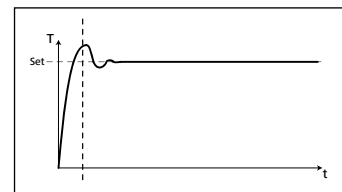
The "Speed limit" menu allows the user to set the desired maximum upper speed limit for the stirrer. The initial setting is the maximum permissible speed of the stirrer. If the user changes this setting, the **Wireless Controller** saves the new value for future stirring tasks. If the "Speed limit" has been changed, then the speed can be adjusted only within the new range.

Control method

The user is able to set the device to three different methods depending on their specific needs. The selected temperature control method is displayed alongside the symbol (11).

**fPID**

**Fast PID:** combines the advantages of the PID and two-point methods (very fast heating and high accuracy; possible overheating).


**aPID**

**Accurate PID:** For use with very sensitive materials which need to be handled carefully.

- Advantages:
- very accurate warming of the liquid
  - no exceeding the target temperature

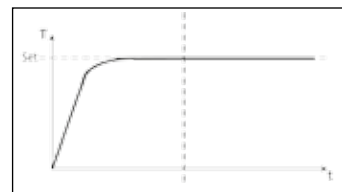
- Disadvantages:
- slow heating operating

**2pt**

**Two-point:** For media which is not too sensitive.

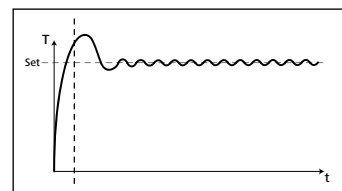- Advantages:
- Heating much more rapid than PID
- Disadvantages:
- Possible overheating of up to 10 K

Limits

In order to ensure that neither the medium nor the carrier fluid overheat, the user can set separate temperature limits.

**Medium limit:**

The "Medium limit" menu allows the user to set the desired maximum temperature limit for the medium.

**Carrier limit:**

The "Carrier limit" menu allows the user to set the desired maximum temperature limit for the carrier fluid.

**Limit information:**

The "Limit information" menu shows the set medium limit and carrier limit.

## Temperature sensor

### **Calibration:**

In order to use tolerances to limit temperature deviations, the user can calibrate the temperature sensor in conjunction with the device.

2-point calibration:

Calibration using two temperatures

3-point calibration:

Most accurate way of calibrating the temperature sensor at three different temperatures

### **Note:**

A calibration carried out by the user is indicated on the display with the symbol **(11)** shown in **red**. The calibration of the PT100 and PT1000.50 double sensor are carried out separately.

### **Reset calibration:**

The menu allows the user to reset the calibration for the temperature sensor.

## Weighing

### Measurement

The weighing module is a function which allows the user to perform simple weighing tasks.

To start the weighing process, press the (G) button on the front membrane or alternatively select "Weighing" from the menu. Press the rotary knob (D) to start the weighing function.

### **Note:**

**The heating and stirring functions must be deactivated.**

It may take several moments for the system to stabilise.

Once the system is stable, a "g" will be shown alongside the measured value.

### Calibration

Place the calibration weight on the mounting plate and wait until the device indicates that the calibration process has been completed.

This process needs to be repeated two or three times before closing calibration mode.

Once the calibration process has been completed successfully, the weighing module is ready for use.

Regularly re-calibrate the device.

## Timer

### Set

The Timer function allows the user to stop the heating process after a certain time has lapsed. The Timer function can be set for up to 99 hours, 59 minutes and 59 seconds.

### **Note:**

**The time starts to count down when starting both functions (heating and/or stirring).**

Once the heating function is activated, the timer function only refers to heating. The stirring function continues to operate as previously, thereby avoiding a delay in boiling in the liquid. The motor is only switched off when the heating function is not active. The stirring function is stopped if the heating function is not started during the timer cycle and the time runs out.

### **Note:**

By entering a value of **00:00:00** (hh:mm:ss), the counter is activated. The counter shows the time elapsed during the current experiment.

### Display

The menu allows the user to specify that the "Timer" is shown on the display. A tick shows that the option is activated.

## pH probe

### Calibration

The device has an integrated pH measuring feature for determining the pH of different liquids.

The pH sensor must be calibrated before being used to attempt a pH measurement.

The calibration is used to adjust the pH probe and the device so that they work together correctly. As part of the process, the neutral and pH gradient are specified for the measurement chain. To complete the calibration, use pH4, pH5 and pH7 buffer solutions in accordance with DIN 19266.

### **Note:**

a pH calibration can only be carried out with an inserted temperature sensor.

- Scroll down to the "pH probe" icon in the menu
- Open the submenu by pressing the rotary knob (D)
- Open the "Calibration" submenu and confirm by pressing the rotary knob (D)
- Place the pH probe and temperature sensor in the required buffer solution.
- Once the value is stable, confirm it by pressing the rotary knob (D).
- Follow the instructions on the display and place the pH probe and temperature sensor in the next buffer solution.
- Confirm once more the value displayed.
- In the case of a 3-point calibration, a third buffer solution must be used.  
pH measurements can be carried out once the calibration has been completed successfully.

### **Note:**

The use of the temperature sensor is limited by the maximum permissible temperature.

Ensure that the temperature of the liquid does not exceed the maximum permissible temperature.

### Display

The menu allows the user to specify that the "pH probe" is shown on the display. A tick shows that the option is activated.

## Mode

The set mode is shown permanently on the display (10).

### Mode A

All the settings are retained when the device is switched off or disconnected from the mains. When the device is switched on, the status of the stirring and heating functions is set to (OFF).

### Mode B

All the settings are retained when the device is switched off or disconnected from the mains. When the device is switched on, the status of the stirring and heating functions is set to whatever it was before the device was last switched off (ON or OFF).

## Mode C

In this mode, the device can be used either with or without a temperature sensor.

When the device is switched on using the main switch, the stirring and heating functions are started in the last set status and at the last set values. The set nominal values in operating modes A and B cannot be modified.

## Display

In this menu the user can specify what information (Torque trend measurement, pH value, Timer or Graph) will be displayed on the main screen. A tick shows that the option is activated.

## Graph

### Axis Assignment

In this menu the user can select 2 options from the 4 options (Temperature, Speed, Torque or pH) for axis assignment. A tick shows that the option is activated.

### Axis Scaling

For axis scaling, the user can choose either the automatic method or the manual method.

## Programs

Under menu "Programs", 5 user-defined speed (rpm)-time profiles can be created. In addition, the user could define the intermittent mode is activated or not in the programs.

If the intermittent mode is activated, the run time/stop time value from "Intermittent Mode" setting will be taken.

**Note:** If user need to activate the intermittent mode in one segment of a program, he should set the "Run time/Stop time" in menu option "Interval", meanwhile activate the "Run/Stop" function in menu option "Intermittent Mode" (see section "Stirring").

Once a program has been selected, the following options are available:

**Start:** Start the selected program upon request.

When the selected program is started by pressing on menu option "Start" with rotary/push knob, the screen allows the user to choose "Infinite loop" or "Loop count". Press "OK" with rotary/push knob to confirm "Infinite loop" and start the program control, or input numbers and then press "OK" with rotary/push knob to confirm "Loop count" and start the program control.

**Edit:** Edit the selected program parameters.

Start to edit the selected program parameters by pressing on menu option "Edit" with rotary/push knob. The user can edit, delete, insert or save one selected program segment in the program.

When user edits the program time for at least one segment, a tick (✓) for respective program will appear.

**Delete:** Delete the selected program.

If a selected program is deleted by pressing on menu option "Delete" with rotary/push knob, all the program parameters will be emptied. The tick (✓) disappears.

**View:** Show details of the selected program.

## Safety

### Time out

The "Time out" function is used to define how long an interruption in communications between the magnetic stirrer and the wireless controller box can be before there is a reaction in the magnetic stirrer.

If the set time limit is exceeded, the pre-selected temperature and speed values become active.

**Note:** The initial setting for the time limit is 30 seconds. The user can set a value of up to 60 minutes for this time limit.

In the "Speed" option, the user can specify a speed that is appropriate and safe for the stirring task. The safe speed applies if there is a communication failure between the stirrer and the **Wireless Controller** or the communication range is exceeded.

**Note:** The initial setting of the safe speed is 50 rpm, and is implemented after expiry of the time limit. The user can set a value of up to 1700 rpm for this speed limit.

If the intermittent mode is activated, the device immediately switches to run at the set safe speed, or the set speed if that is less than the safe speed.

In the "Temperature" option, the user can specify a temperature that is appropriate and safe for the heating task. The safe temperature applies if there is a communication failure between the stirrer and the **Wireless Controller** or the communication range is exceeded.

**Note:** The setting range for the safe temperature is 0 - 340 °C.

### Password

In the "Password" menu, the user can protect the **Wireless Controller** settings using a password (factory setting: 000).

### Safe temp confirmation

This option requires the user to accept the hot plate safe temperature value during the start-up.

### Er5

The time is taken to recognise that the sensor has not been placed in the medium when the heating is switched on.

**Note:** The user can set a value from 1 to 30 min for this time limit depending on the application.

If time limit is set to 0 min, the Error 5 detection is disabled.

This function will only be active if:

- Sensor temperature is < 50 °C
- Difference target temperature / sensor temperature > 5 K

## Settings

### Languages:

The "**Languages**" option allows the user to select the desired language by turning and pressing the rotary/push knob (D). A tick indicates the language that is set for the system.

### Units:

The "**Units**" option allows the user to select the desired unit for displaying the temperature. The choice is between "**°C**" or "**°F**" and is made by turning and pressing the rotary/push knob (D). A tick indicates the unit that is set for the system.

### Display:

The "**Display**" option allows the user to change the background color between "Black" and "White", and display "Firmware Update Information". A tick shows that the option is activated.

### Sound:

The "**Sound**" option allows the user to activate/deactivate the key tone. A tick shows that the option is activated.

### Factory Settings:

Select the "**Factory Settings**" option by turning and pressing the rotary/push knob. The system requests confirmation to restore the factory settings. Pressing the "OK" button resets all the system settings to the original standard values set at dispatch from the factory (see "Menu structure").

### Communication:

#### Bluetooth

The "**Bluetooth**" option allows the user to activate/deactivate the "**Bluetooth**" function. A tick shows that the option is activated.

The device allows the user to establish a connection to a wireless controller box.

The device can then be operated via a wireless controller box within a range of 10 m (without obstructions).

Please refer to the separate instruction manual for the wireless controller box for more details.

#### Eurostar control

The new safety feature of the RET monitors the behavior of a connected Eurostar and stops the heating process when the Eurostar stirring speed falls below a specific minimum speed. The RET also stops the heating process whenever the Eurostar is in an error state or the cable connection between the RET and the Eurostar gets lost.

The Eurostar can be connected with the RET control either via RS232 or USB. The used interface has to be selected in the menu.

**Main Menu → Settings → Communication → Eurostar control**  
**Important information when the USB interface is used for the communication between the RET and the Eurostar:**

The USB cable has always be connected with the RET first and then with the Eurostar.

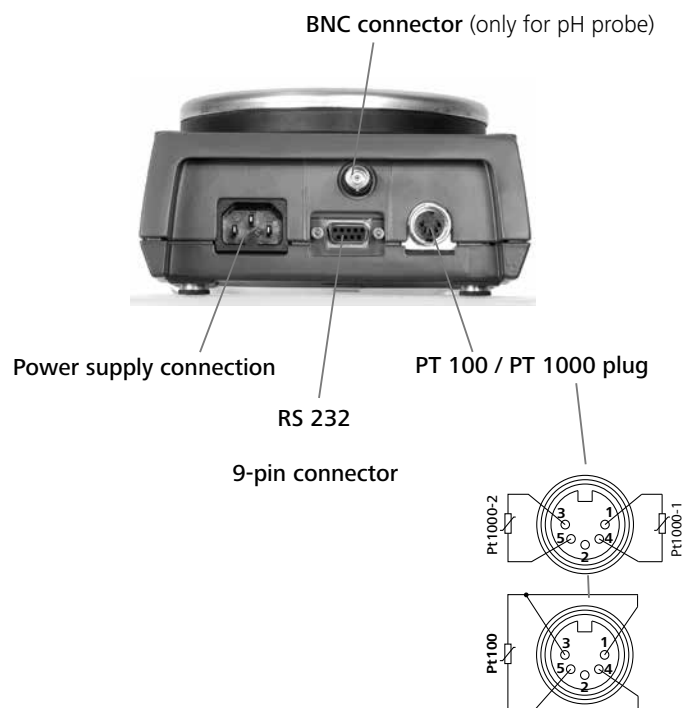
For the stirrer speed of the Eurostar has to be a minimum speed set. When the actual value of the stirrer speed falls below the minimum limit the heating function of the RET will be stopped and the RET will be set in an error state.

The new menus are only in English.

#### Information:

The "**Information**" option offers the user an overview of the most important system settings of the stirrer RET control-visc.

## Connections and interfaces

## Serial RS 232 interface and USB

### USB interface

The Universal Serial Bus (USB) is a serial bus for connecting the device to the PC. Equipped with USB devices can be connected to a PC during operation (hot plugging). Connected devices and their properties are automatically recognized. Use the USB interface in conjunction with labworldsoft® for operation in "Remote" mode and also to update the firmware.

### Installation

First, download the latest driver for IKA® devices with USB interface from <http://www.ika.com/ika/lws/download/usb-driver.zip> and install the driver by running the setup file. Then connect the IKA® device through the USB data cable to the PC. The data communication is via a virtual COM port. Configuration, command syntax and commands of the virtual COM ports are as described in RS 232 interface.

### Configuration

The RS 232 serial interface can be used to operate the device using a computer and a suitable user program (labworldsoft).

**In order to increase safety when operating the RET using a PC, a watchdog function can be activated which monitors the continuous data flow (see section entitled: "Watchdog function").**

- The functions of the interface circuit between the laboratory device and the automation system are a selection from the signals specified in the EIA standard RS232 as per DIN 66020 Part 1. The assignment of the different signals can be seen in the image.
- Standard RS 232, corresponding to DIN 66259 Part 1 is valid for the electronic characteristics of the interface circuits and assignment of signal states.
- Transmission process: asynchronous character transmission in start-stop operation.
- Transmission type: full duplex
- Character format: character composition according to data format in DIN 66022 for start-stop operation. 1 start bit, 7 character bits, 1 parity bit (even), 1 stop bit.
- Transmission speed: 9,600 bit/s
- Data flow control: hardware handshake RTS/CTS
  - RTS: (PIN 7) LOW (positive voltage)  
/ COMPUTER can transmit
  - RTS: (PIN 7) HIGH (negative voltage)  
/ COMPUTER cannot transmit

CTS: (PIN 8) LOW (positive voltage)  
/ COMPUTER ready-to-receive  
CTS: (PIN 8) HIGH (negative voltage)  
/ COMPUTER not ready-to-receive

- Access method: data transmission from the device to the computer only occurs after a request from the computer.

### Command syntax

The following points should be noted for the instruction set:

- Commands are generally sent from the computer (master) to the lab device (slave).
- The lab device only responds to requests from the computer.
- Commands and parameters, as well as consecutive parameters, must be separated by at least one space (code: hex 0x20).
- Each individual command including parameters and data and all responses are completed with CR LF (code: hex 0x0D and 0x0A) and can have a maximum length of 10 characters.
- The dot is used for decimal separators in a floating-point value (code: hex 0x2E).

The details given above generally comply with the recommendations of NAMUR (NAMUR recommendations for the design of electrical plug-in connectors for analogue and digital signals in laboratory MSR devices. Rev. 1.1).

### Summary of available NAMUR commands

Abbreviations used:

- X,y= numbering parameter (integer)
- m = variable value, integer
- n = variable value, floating-point number
- X = 1 medium temperature(external temperature sensor)
- X = 2 hot plate temperature
- X = 3 hot plate safety temperature
- X = 4 speed
- X = 5 viscosity trend
- X = 7 heat transfer medium temperature
- X = 80 pH value
- X = 90 Weight value
- X = 54 Error 5 response time in seconds (180 ≤ n ≤ 1200)
- X = 55 Intermittent mode cycle time in seconds  
(10 ≤ n ≤ 600)
- X = 56 Intermittent mode, duration of interruption in seconds  
(5 ≤ n ≤ 60)

NAMUR commands	Function	Display (additional)
IN_NAME	Title request	
IN_PV_X X=1;2;3;4;5;7;80;90	Current value reading	
IN_SOFTWARE	Software ID number, date, version request	
IN_SP_X X=1;2;3;4;7;8;12;42;54;55;56;	Set target value reading	
IN_TYPE	Lab device identification request	
OUT_NAME name	Output of identification name. (Max. 6 characters; default: IKARET)	
OUT_SP_12@n	Setting WD safety limit temperature with set value echo	
OUT_SP_42@n	Setting WD safety limit speed with set value echo	
OUT_SP_X n X=1;2;4;7;54;55;56	Setting of target value to n	

OUT_WD1@m	Watchdog mode 1: if event WD1 should occur, the heating and stirring functions are switched off and ER 2 is displayed. Set watchdog time to m (20 - 1,500) seconds, with watchdog time echo. This command launches the watchdog function and must be transmitted within the set watchdog time.	
OUT_WD2@m	Watchdog mode 2: if event WD2 should occur, the speed target value is changed to the WD safety speed limit and the temperature target value is changed to the WD safety temperature limit value. The warning WD is displayed. The WD2 event can be reset with OUT_WD2@0 - this also stops the watchdog function. Set watchdog time to m (20 - 1,500) seconds, with watchdog time echo. This command launches the watchdog function and must be transmitted within the set watchdog time.	
RESET	Switches off the device function.	
START_X X=1;2;4;5;7 X=80; 90	Switching on of device - (remote) function Activation of the previous watchdog function. If no target value is entered within 60 seconds, the (old function) heating is switched off.	Remote
STOP_X X=1;2;4;5;7 X=80; 90	Switching off of device - (remote) function The variables set with OUT_SP_X remain saved. Includes command RMP_STOP.	Remote

### “Watchdog” functions; monitoring of the serial data flow

If, once this function has been activated (see NAMUR commands), there is no retransmission of the command from the computer within the set time (“watchdog time”), the heating and stirring functions are switched off in accordance with the set “watchdog” function or are changed to the set target values.

The data transmission may be interrupted by, for example, a crash in the operating system, a power failure in the PC or an issue with the connection table between the computer and RET control-visc.

#### “Watchdog” - mode 1

If there is an interruption in data communications (longer than the set watchdog time), the heating and stirring functions are switched off and ER 2 is displayed.

#### “Watchdog” - mode 2

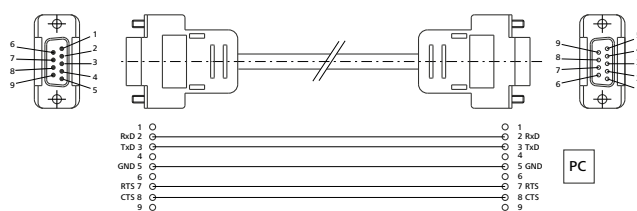
If there is an interruption in data communications (longer than the set watchdog time), the speed target value is changed to the WD safety speed limit and the temperature target value is changed to the WD safety temperature limit value. The warning WD is displayed.

### Connection possibilities between the RET control-visc and external devices

The following adapters and adapter cables are available from IKA® for connecting the RET control-visc to external devices.

#### PC 1.1 cable

The PC 1.1 cable is used to connect the 9-pin plug to a computer.

## Accessories

- Magnetic stirring bar:
  - ø 6 mm; max. length 15 mm
  - ø 7 mm; max. length 60 mm
  - ø 10 mm; max. length 80 mm
- RS 1 Set of magnetic stirring bars
- RSE Stirring bar remover
- H 1000 Beaker 1 l
- H 1500 Beaker 1.5 l
- H 3000 Beaker 3 l
- H 5000 Beaker 5 l
- H 8000 Beaker 8 l
- H 135.3 Flask carrier
- H 135.310 Quarter 20 ml reaction vessel
- H 135.311 Quarter 30 ml reaction vessel
- H 135.312 Quarter 40 ml reaction vessel
- H 135.313 Quarter 4 ml reaction vessel
- H 135.314 Quarter 8 ml reaction vessel
- H 135.315 Quarter 16 ml reaction vessel
- H 135.4 Reaction block for 100 ml round flask
- H 135.410 Insert for 10 ml round flask
- H 135.411 Insert for 25 ml round flask
- H 135.412 Insert for 50 ml round flask
- H 135.5 Reaction block for 500 ml round flask
- H 135.510 Insert for 200 ml round flask
- H 135.511 Insert for 250 ml round flask
- H 135.512 Insert for 300 ml round flask
- H 135.6 Reaction block for 1,000 ml round flask
- H 135.610 Insert for 500 ml round flask
- H 16 V Support rod
- H 16.3 Extension
- H 38 Holding rod
- H 44 Cross sleeve
- PT 100.50 Temperature sensor
- PT 100.51 Temperature sensor, glass coated
- PT 1000.50 Double temperature sensor
- PT 1000.51 Double temperature sensor, glass coated
- RET control-visc WiCo

for further accessories see [www.ika.com](http://www.ika.com).

## Troubleshooting

The device will indicate that there is a fault by displaying an error message "Er X" (see table) on the screen.  
Heating will be switched off permanently (exception: WD - see table).

The stirring function continues to operate at the pre-set speed if there is a fault in the heating circuit (exception: WD - see table).  
In order to reset the device after a fault, it may need to be disconnected from the power supply for several seconds. The period the device is switched off may also need to be longer if the temperatures need to be counterbalanced (residual heat display may no longer be active; see "Residual heat display").

The safety circuit test takes place within 20 seconds of switching on the device; after the test, the other monitoring functions are activated.

**If a fault cannot be eliminated directly, you must perform a RESET** (see "Settings" menu). **If the faults still cannot be eliminated, the appliance must be inspected by a technical service technician.**

### Error messages and troubleshooting table

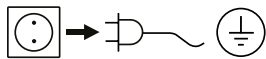
Fault	Description	Causes	Effect	Solutions
Er 2	In remote operation (PC) mode with activated watchdog function in mode 1: no communication between PC and RET control-visc	<ul style="list-style-type: none"> <li>PC does not transmit any data within the set watchdog time</li> <li>Connection to PC interrupted</li> </ul>	Heating switched off Motor switched off	<ul style="list-style-type: none"> <li>Change watchdog time</li> <li>Transmit data from PC within set watchdog time (OUT_WDx@m)</li> <li>Check cable and plug</li> </ul>
Er 3	Device internal temperature exceeds 80 °C	<ul style="list-style-type: none"> <li>Heat accumulation between heating plate and housing</li> <li>Permitted ambient temperature exceeded</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Switch device off, leave to cool and switch on again</li> <li>Change experiment</li> <li>Observe maximum permissible ambient temperature</li> </ul>
Er 4	Difference between target and actual speed exceeds 300 rpm	<ul style="list-style-type: none"> <li>Motor blocked or overloaded</li> </ul>	Heating switched off Motor switched off	<ul style="list-style-type: none"> <li>Reduce load torque or use smaller magnetic rods</li> <li>Reduce target speed</li> </ul>
Er 5	No temperature increase in sensor at continuous temperature difference (selected time in menu)	<ul style="list-style-type: none"> <li>Sensor not in medium</li> <li>Volume of medium to be measured too large</li> <li>Heat conductivity of medium to be measured too low</li> <li>Heat conductivity of the vessel is too low</li> <li>In the case of indirect heating, the overall heat conductivity resistance is too large</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Place the sensor in the medium</li> <li>Reduce the volume of the media</li> <li>Use a carrier fluid with better heat conductivity properties</li> <li>Replace the glass vessel with a metal pot</li> <li>Increase the "Time-out" period</li> </ul>
Er 6	Interruption in safety circuit	<ul style="list-style-type: none"> <li>Defective connection cable</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Replace cable</li> </ul>
Er 13	SAFE TEMP P adjustable safety circuit defective	<ul style="list-style-type: none"> <li>Target/current different of the adjustable safety circuit for minimum temperature monitoring</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>After switching on, change the SAFE TEMP to a different value; if this solves the issue, the previous value can be reset by switching the device off and on again</li> </ul>
Er 14	Short circuit in temperature sensor 1	<ul style="list-style-type: none"> <li>Short circuit in temperature sensor plug</li> <li>Short circuit in the cable or temperature sensor</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Check the plug</li> <li>Replace the temperature sensor</li> </ul>
Er 21	Fault during heating plate safety test	<ul style="list-style-type: none"> <li>Safety relay does not open</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Switch device off, leave to cool and switch on again</li> </ul>
Er 22	Fault during heating plate safety test	<ul style="list-style-type: none"> <li>S_CHECK cannot generate H_S_TEMP</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Switch device off, leave to cool and switch on again</li> </ul>
Er 24	The temperature of the heating plate exceeds SAFE TEMP H (Hotplate)	<ul style="list-style-type: none"> <li>SAFE TEMP H (Probe) has been set to lower than the current temperature of the heating plate</li> <li>Disconnection of heating plate control temperature sensor</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>Leave the heating plate to cool</li> <li>Set the SAFE TEMP H (Hotplate) higher</li> </ul>

Fault	Description	Causes	Effect	Solutions
Er 25	Heating switching element monitoring	<ul style="list-style-type: none"> <li>• Heater control circuit switch (TRIAC) short-circuited</li> <li>• Safety relay has interrupted the heating circuit</li> <li>• Heater or the supply line is disconnected</li> <li>• Interruption in heating plate safety temperature sensor</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>• Switch device off, leave to cool and switch on again</li> </ul>
Er 26	Difference between internal heating plate control and safety temperature sensor too large	<ul style="list-style-type: none"> <li>• Irregular temperature distribution across heating plate due to sporadic heat dissipation</li> <li>• Defective control or safety temperature sensor</li> </ul>	Heating switched off	<ul style="list-style-type: none"> <li>• Switch device off, leave to cool and switch on again</li> <li>• Ensure regular heat dissipation when using metal blocks, etc. on the flat surface of the heating plate</li> </ul>
Er 31	Defective monitoring of the heating control circuit (TRIAC)		Heating switched off	<ul style="list-style-type: none"> <li>• Contact customer service</li> </ul>
Er 41	Bar decoupling	<ul style="list-style-type: none"> <li>• The bar has broken away four times in a row without an explicit change in the target value</li> </ul>	Target speed reduced by at least 300 rpm Heating switched off	<ul style="list-style-type: none"> <li>• Select a lower target speed</li> <li>• Select a smaller stirring bar</li> </ul>

## Maintenance and cleaning

The device should operate without the need for maintenance, however it is subject to natural wear and tear on parts and their statistical failure rate.

### Cleaning

Remove the device from the mains before cleaning.

Only use cleaning materials recommended by IKA®.

Dirt	Cleaning agent
Dyes	Isopropyl alcohol
Building materials	Water containing detergent/isopropyl alcohol
Cosmetics	Water containing detergent/isopropyl alcohol
Food	Water containing detergent
Fuels	Water containing detergent

- Ensure no liquid enters the device during cleaning.
- Wear protective gloves when cleaning the device.
- Please consult with IKA® before using any cleaning or decontamination methods not specifically recommended.

### Ordering spare parts

When ordering spare parts, please make sure to indicate the following:

- device type
- device manufacturing number; see rating plate
- software version (second value displayed when switching on the device)
- position number and description of spare part; see [www.ika.com](http://www.ika.com).

### Repairs

**Please only send devices in for repair that have been cleaned and are free of materials which might present health hazards.**

For this, use the “**certificate of compliance**” form which you can obtain from IKA® or can download a version for printing from the IKA® website at [www.ika.com](http://www.ika.com).

If your appliance requires repair, return it in its original packaging. Storage packaging is not sufficient when sending the device - also use appropriate transport packaging.

## Technical data

Rated voltage	or or	VAC VAC VAC	220...230 ± 10% 115 ± 10% 100 ± 10%
Frequency		Hz	50 / 60
Power input		W	650
Display			TFT
Timer		h	99 h 59 min 59 sec
Stirring			
Motor:			ball bearing mounted, brushless EC motor

Number of stirring positions		1
Max. motor power consumption	W	22
Max. motor output	W	12
Speed range	rpm	0; 50 to 1700
Target speed setting resolution	rpm	1
Actual speed display resolution	rpm	1
Speed variation (without load, 115/230 V, at 1700 rpm, RT 25 °C)	%	2
Stirring capacity (H <sub>2</sub> O)	l	20
<u>Heating</u>		
Self-heating of the hotplate (RT:22 °C/duration:1 h)	°C	35
Heating plate material		Stainless steel 1.4301
Heating plate diameter	mm	135
Heat output (100/115/230 V)	W	600 ± 5%
Heating rate (1 l H <sub>2</sub> O in H15)	K/min	7
Temperature range	°C	RT ... 340
Temperature medium max. (dep. on vessel)	°C	265
Target temperature setting resolution	K	0.1
Actual temperature display resolution - of the medium temperature - of the carrier fluid temperature	K K	0.1 1
Adjustable heating plate safety circuit	°C	50 - 380 ( ± 5 °C)
Safety circuit setting resolution	K	1
Temperature sensor in medium		PT 100 or 2x PT 1000, DIN EN 60751 Class A
Temperature sensor deviation (PT 100 or 2x PT 1000, DIN EN 60751 Class A)	K	<= ± [0.15 + 0.002 °(T)]
Deviation in section of temperature measurements	K K	<= ± 0.3 < 100 °C <= ± 0.5 > 100 °C
Deviation in section of temperature measurements + sensor	K K	<= ± 0.45 at 0 °C <= ± 0.85 at 100 °C
Deviation in heating plate temperature	K	± 3 without vessel; heating plate at 100 °C
Control hysteresis with temperature sensor (500 ml water, 600 ml glass beaker, 30 mm bar, 800 rpm, PT 100.70, 50 °C)	K	± 0.2
Heating plate control hysteresis	K	± 3 without vessel; heating plate at 100 °C
Permissible ambient temperature	°C	+ 5 to + 40
Permissible relative humidity	%	80
Permissible operating time	%	100
Protection class according to DIN EN 60529		IP 42
Weighing range	g	10 - 5000
Load capacity weighing function max.	g	5000
Weighing range tolerance	g	< 500: ± 1 > 500: ± 5
Level of contamination		2
Overvoltage category		II
Protection class		1 (protective earth)
Use above sea level	m	max. 2000
Dimensions (W x D x H)	mm	160 x 270 x 85
Weight	kg	2.7

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**IKA® - Werke GmbH & Co.KG**