

# dataLink™

Traceability system for easySpiral® platers  
and Scan® automatic colony counters

## 100% traceability

from plating to colony counting



# interscience



easy**Spiral** Dilute®  
serial dilutor and automatic plater



**Scan**®  
automatic colony counter

**dataLink™** is a full traceability system for the microbiological analysis of a sample, from plating to colony counting with **easySpiral®** automatic platers and **Scan®** automatic colony counters.

All plating data (volume, dilution, time, mode ...) is stored on a Datamatrix or QR Code label placed on the Petri dish. No database is needed.

After incubation, the Datamatrix label is scanned and read by the **Scan®** software and retrieves all the sample information, from the plating to the colony counter settings.

Simply press "COUNT". There is no more time-consuming data double input and potential errors. Export the results, pictures and comments to LIMS\*, Excel™, PDF, JPEG.

### ► **Secured analyses**

- Full traceability
- Error-free settings

### ► **Time-saving**

- No manual data input
- Automatic settings of the Scan® colony counter

### ► **User-friendly**

- Stand-alone system
- No database or network use

## STEP 1


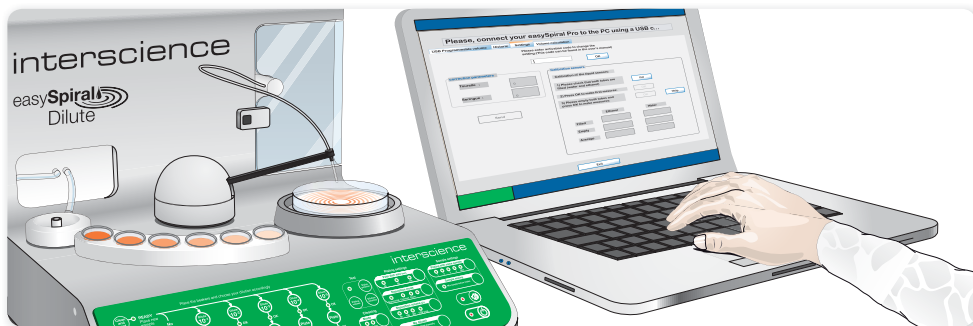


Plate with easy**Spiral**® Pro or Dilute. easy**Spiral**® software collects the plating data.

## STEP 2

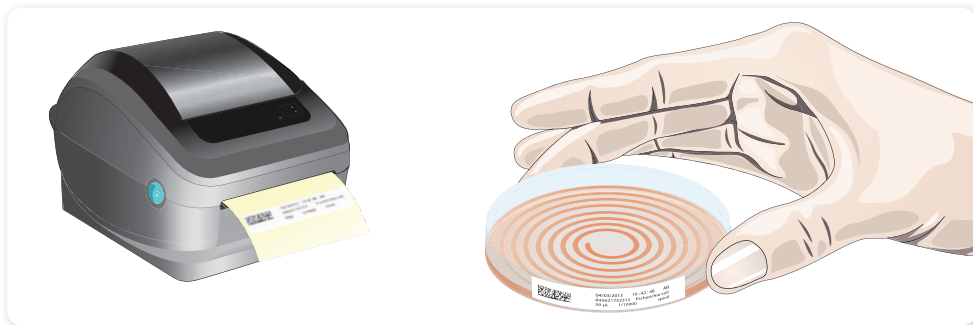

Print the label with Datamatrix code. Stick the label on the plated Petri dish and place in the incubator.

■■■ Incubation 24-72 h

## STEP 3



Once the colonies have grown, scan the Datamatrix code. The **Scan**® colony counter automatically adjusts its settings thanks to the Datamatrix label's data. Click on "COUNT". Export the data.

\* Please check LIMS compatibility

# 5 label models:




## Bar-code 39




## Bar-code 128



## Datamatrix

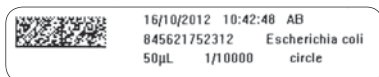

## QR Code

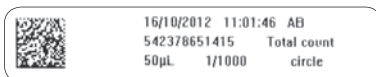

## Number




Label dimensions: 45 x 10 mm

## dataLink™ pack:



Ref. 410 100

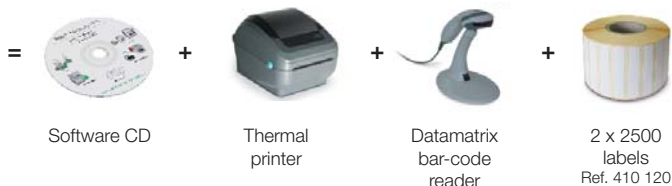

Your local distributor

## interscience INTERNATIONAL

30, ch. Bois Arpents - 78860 St Nom - FRANCE  
T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03  
Email: [info@interscience.com](mailto:info@interscience.com)

## interscience USA & CANADA

32 Cummings Park - Woburn, MA 01801 - USA  
T: +1 781 937 0007 - F: +1 781 937 0017  
Email: [sales.usa@interscience.com](mailto:sales.usa@interscience.com)

## interscience CHINA

上海徐汇区吴兴路277号锦都大厦718室-200030  
电话: +86(0)21-64739390 - 手机: +86 189 3097 0733  
邮址: [sales.china@interscience.com](mailto:sales.china@interscience.com)

## interscience SOUTH-EAST ASIA

541 Orchard Rd -09-01 Liat Towers - SINGAPORE 238881  
T: +65 6933 1389 - M: +65 8549 1217  
Email: [sales.asia@interscience.com](mailto:sales.asia@interscience.com)

[www.interscience.com](http://www.interscience.com)