
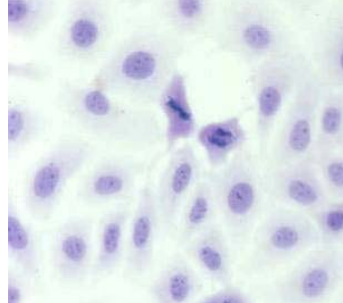


Single Cell Isolation Using Glass Chips

Protocol

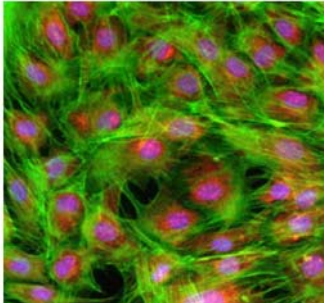

CORNING

*John A. Ryan, Ph.D.
Corning Incorporated
Life Sciences
900 Chelmsford St
Lowell, MA 01851*

Introduction

Glass chips have two advantages over other clonal isolation techniques. First, visual observation is generally easier; the presence of a single cell can be easily verified. Second, if you have a steady hand, a fairly large number of chips can be picked in a short time period. A disadvantage of the method is that when working with primary cultures or cell lines that have a low cloning efficiency, many chips must be picked before one will be found that will develop into a colony. Another disadvantage is that preparing the chips is tedious.

Materials

Supplies

1. Watchmaker forceps
2. Stainless steel U.S. Standard soil sieves (1 mm and .45 mm diameters)
3. 24 well microplates (Corning[®] Cat. #3526 or 3524)
4. 60 mm culture dishes (Corning Cat. #430166)
5. Glass coverslips (Corning Cat. #2875-25 or similar)
6. Mortar and pestle (must be very clean)
7. Concentrated nitric acid
8. Glass beaker (1 L)
9. Glass tubes with metal caps

Medium and Solutions

1. Calcium-and magnesium-free phosphate buffered saline (CMF-PBS) for dissociating cultures

2. 0.25% Trypsin in CMF-PBS (or other enzyme solution suitable for cells being picked)
3. Growth medium appropriate for cells being picked.
4. Cell culture for harvesting

Procedure

1. Break coverslips into 1 mm² to 2 mm² fragments. (This can be done by a variety of methods. One suggestion is using a mortar and pestle.)
2. The fragments are then sifted through two stainless steel U.S. Standard soil sieves (1 mm and .45 mm diameters). The fragments retained on the .45 mm screen are then washed. The larger fragments can be collected and broken into smaller fragments and rescreened.

SAFETY PRECAUTION: THE FOLLOWING STEP MUST BE DONE UNDER A FUME HOOD.

3. Put the coverslip fragments into a large PYREX[®] beaker and add equal volumes of water and nitric acid (add acid to water). Then bring to a fuming point (60 to 70°C) on an electric hot plate in a fume hood for 30 minutes. Remove beaker from hot plate to cool. When cool, decant the acid into the hood's cup sink with water running; rinse the chips thoroughly under cold running water for no less than three hours, then rinse six times in distilled water.
4. Transfer the sized coverslip fragments to glass tubes with metal closures and dry heat sterilize at 180°C for 3 hours.
5. Add 50 to 200 coverslip fragments to each of six 60 mm plastic tissue culture dishes and distribute evenly across bottoms.
6. Carefully add 3 ml of growth medium to each plate. Avoid floating the glass fragments.
7. Prepare cell suspension by the gentlest means possible.
8. Inoculate (using 1 mL of medium) duplicate dishes containing glass fragments with 1, 2, and 5 x 10³ cells per dish. It is important to do a range of cell concentrations. If the initial cell concentration is too high most of the chips will have more than one cell. If it is too low then very few if any chips will have a cell.
9. After cells have had sufficient time to attach, (1 to two hours) look for coverslip fragments with a single adherent cell. Start with the cell concentration that has the best ratio of chips with single cells.
10. Using watchmaker forceps transfer those fragments with a single cell to separate wells of a 24 well plate containing 1 mL medium/well (Figure 1). Use conditioned medium if possible; it will increase the cloning efficiency as well as help the cells grow faster. Make sure the cell is on the top side of the glass chip after it is placed into the well or else the cell may be unable to divide.
11. Once the surface of the glass fragments is overgrown with cells (usually 5-10 days), it is best to disperse the colony by enzymatic treatment. The cells can then be replated in a similar culture vessel. When left alone on the glass

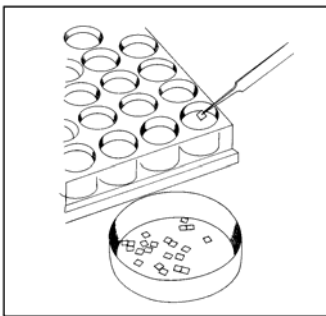


Figure 1. It takes a steady hand to pick up the glass chips and is initially quite challenging. But with practice it is possible to find and remove chips quickly and efficiently.

fragments, colonies will migrate onto the dish surface, but growth is usually slower in this manner.

Acknowledgments

This protocol has evolved from protocols developed for cell culture training courses at the former W. Alton Jones Cell Science Center in Lake Placid, New York; Manhattan College, New York City; and the University of Connecticut, Storrs, Connecticut. I would like to thank all of my colleagues and students who, over the years, have contributed ideas and suggestions to its development.

References

1. Martin, G. M. (1973). **Dilution Plating on Coverslip Fragments**. In: Tissue Culture Methods and Applications. P. F. Kruse and M. K. Patterson Eds. Academic Press, New York, pages 264-266.
2. Freshney, R. I. (2000). **Culture of Animal Cells: A Manual of Basic Technique**. Fourth edition, Wiley-Liss, Inc. New York, pages 194 – 213.
3. Reid, L. (1979). **Cloning**. In: Methods in Enzymology, Vol. 58: Cell Culture. W. B. Jakoby and I. H. Pastan, Eds. Academic Press, New York, pages 152 – 164.

Please visit the Corning web site (www.corning.com/lifesciences) to download these and other useful cell culture articles and protocols

[Mycoplasma Detection using DNA Staining](#)

This protocol provides a detailed procedure for using Hoechst stain #33258 to stain the DNA in cell monolayers for detecting mycoplasmas or other prokaryotic organisms.

[Subculturing Monolayer Cell Cultures](#)

This protocol provides a detailed procedure for subculturing attached cells using enzymatic dissociation. It explains the theory behind some of the key steps in the process.

[Cryogenic Preservation and Storage of Animal Cells](#)

This protocol provides a detailed procedure for cryogenically storing a variety of animal cells.

[Calcium Phosphate Transfections in Corning® CellSTACK® Culture Chambers](#)

This protocol is a generalized procedure for calcium phosphate based transfection in Corning CellSTACK Chambers. It is intended as a starting point for individuals wishing to transfect large numbers of cells in the Corning CellSTACK chamber and can be adapted for each size.

[Single Cell Cloning by Serial Dilution](#)

This protocol describes a simple approach using a 96 well plate to isolate colonies derived from single cells.

Clonal Growth of Cells in Semisolid Media

This protocol provides a detailed procedure for growing animal cells in agar-based media for either transformation assays or cloning.

Use of Corning Cloning Cylinders for Harvesting Cell Colonies

This protocol describes how to isolate and develop a population of cells (a clone) that are the descendants of a single cell.

Harvesting Cells from CellSTACK® Chambers

This protocol gives instructions for harvesting cells from 10-layer CellSTACK chambers and discusses some of the options and products that can make this process easier.

For additional product or technical information, please e-mail us at CLStechserv@corning.com, visit our web site www.corning.com/lifesciences or call 1-800-492-1110. Outside the United States call 978-635-2200.

CORNING

Corning Incorporated *Life Sciences*

Tower 2, 4th Floor
900 Chelmsford St.
Lowell, MA 01851
t 800.492.1110
t 978.442.2200
f 978.442.2476

www.corning.com/lifesciences

Worldwide Support Offices

ASIA/PACIFIC

Australia
t 61 2-9416-0492
f 61 2-9416-0493

China
t 86 21-3222-4666
f 86 21-6288-1575

Hong Kong
t 852-2807-2723
f 852-2807-2152

India
t 91 11 341 3440
f 91 11 341 1520

Japan
t 81 (0) 3-3586 1996/1997
f 81 (0) 3-3586 1291/1292

Korea
t 82 2-796-9500
f 82 2-796-9300

Singapore
t 65 6733-6511
f 65 6735-2913

Taiwan
t 886 2-2716-0338
f 886 2-2716-0339

EUROPE

France
t 0800 916 882
f 0800 918 636

Germany
t 0800 101 1153
f 0800 101 2427

United Kingdom
t 0800 376 8660
f 0800 279 1117

The Netherlands
t 31 (0) 20 659 60 51
f 31 (0) 20 659 76 73

All other European Countries
t 31(0) 20 659 60 51
f 31(0) 20 659 76 73

**LATIN AMERICA
Brasil**
t (55-11) 3089-7400
f (55-11) 6845-2236

Mexico
t (52-81) 8313-8400
f (52-81) 8313-8589

Corning, PYREX and CellSTACK are registered trademarks of Corning Incorporated, Corning, New York.

Corning Incorporated, One Riverfront Plaza, Corning, NY, 14831-0001